

# Package ‘AffyCompatible’

October 7, 2014

**Type** Package

**Title** Affymetrix GeneChip software compatibility

**Version** 1.24.1

**Author** Martin Morgan, Robert Gentleman

**Maintainer** Martin Morgan <mtmorgan@fhcrc.org>

**Depends** R (>= 2.7.0), XML (>= 2.8-1), RCurl (>= 0.8-1), methods

**Imports** Biostrings

**Description** This package provides an interface to Affymetrix chip annotation and sample attribute files. The package allows an easy way for users to download and manage local data bases of Affymetrix NetAffx annotation files. The package also provides access to GeneChip Operating System (GCOS) and GeneChip Command Console (AGCC)-compatible sample annotation files.

**License** Artistic-2.0

**biocViews** Infrastructure, Microarray, OneChannel

## R topics documented:

|                                  |    |
|----------------------------------|----|
| AffyCompatible-package . . . . . | 2  |
| accessors . . . . .              | 2  |
| AffyCompatible-class . . . . .   | 3  |
| Internal-class . . . . .         | 4  |
| NetAffxResource . . . . .        | 4  |
| NetAffxResource-class . . . . .  | 6  |
| readArr . . . . .                | 8  |
| readMage . . . . .               | 9  |
| readNetAffx . . . . .            | 10 |
| readXml . . . . .                | 11 |
| xclass . . . . .                 | 12 |

|  |    |
|--|----|
| xmlValue.XMLAttributeValue . . . . .     | 13 |
| zzz-ARRCompatibility-class . . . . .     | 13 |
| zzz-DTTCompatibility-class . . . . .     | 14 |
| zzz-NetAffxCompatibility-class . . . . . | 16 |

|              |           |
|--------------|-----------|
| <b>Index</b> | <b>18</b> |
|--------------|-----------|

---

AffyCompatible-package

*Affymetrix chip annotation and sample attribute management*

---

## Description

This package manages Affymetrix NetAffx annotation files, and parses MAGE and ARR sample attribute files into R data structures.

NetAffx provides a way to manage the diverse annotation files provided by Affymetrix for their arrays.

ARR files provide attribute information about samples used in an experiment. These files are created, for instance, by the GeneChip Command Console or compliant applications, or

MAGE files also provide information about samples. Support in this package is for MAGE produced by Affymetrix GCOS compliant applications, and emphasizes sample description and processing.

## Details

Package details are available in the DESCRIPTION file of the package, or with the R command `packageDescription("AffyCompatible")`

Overviews of how to use the functionality of this package are available in the package vignettes. See `vignette(package="AffyCompatible")`.

## Author(s)

Written by Martin Morgan.

Maintainer: Martin Morgan <martin.t.morgan@gmail.com>

---

accessors

*Access NetAffxResource-class content*

---

## Description

These functions exist to access the information represented in objects of `linkS4class{NetAffxResource}` and related classes.

**Usage**

```
annotationsFile(x)
directory(x)
password(x)
user(x)
```

**Arguments**

x                    An object for which the accessor is defined.

**Value**

See, e.g., `help(directory, NetAffxResource-method)` for values returned by specific methods.

**Author(s)**

Martin Morgan

**See Also**

[NetAffxResource](#)

---

AffyCompatible-class    *Virtual classes defining the "AffyCompatibility" hierarchy*

---

**Description**

The classes `AffyCompatible`, `DTTCompatibility`, `ARRCompatibility`, and `NetAffxCompatibility` define the overall class hierarchy in this package. Information about these classes is not usually relevant to the user.

**Objects from the Class**

These are virtual classes: no objects may be created from them.

**Methods**

`show` signature(object = "AffyCompatible"): for each slot, provide the slot name, and then invoke `show` for the slot content.

**Author(s)**

Martin Morgan

**Examples**

```
showClass(getClass("NetAffxCompatibility",
  where=getNamespace("AffyCompatible")))
```

---

Internal-class                    *(Internal) Represent 'typed associations' from XML documents*

---

### Description

This internal class represents 'typed associates' (vectors of identically-typed objects) that appear as vectors to the R user.

### Objects from the Class

Objects are created internally, during XML document parsing.

### Slots

`.Data`: Object of class "list" containing the association objects  
`.extends`: Object of class "character" and length 1 describing the class each element in `.Data` must extend

### Extends

Class "list", from data part. Class "vector", by class "list", distance 2.

### Methods

[ signature(x = ".TypedAssociation", i = "numeric", j = "missing"): Subset the typed association, returning a typed association with elements i.  
 [ signature(x = ".TypedAssociation", i = "ANY", j = "ANY"): This method is present to stop access to subsetting methods for class list; it has no user-visible functionality.  
**show** signature(object = ".TypedAssociation"): Compactly display the typed associate, e.g., for `.extends=Foo` and length 5, display `Foo(5)`.

### Author(s)

Martin Morgan

---

NetAffxResource                    *Create and access a NetAffxResource database*

---

### Description

Use this function to create and access a NetAffxResource database.

**Usage**

```
NetAffxResource(user = character(0),  
  password = character(0),  
  affxLicence = "FHCRC0607",  
  directory = tempdir(), ...)
```

**Arguments**

|             |   |
|-------------|---|
| user        | Length 1 character vector containing the Affymetrix user name authorized to access the NetAffx resource.  |
| password    | Length 1 character vector containing the Affymetrix password of the user.   |
| affxLicence | Length 1 character vector containing the Affymetrix licence (typically application specific, i.e., the specified licence covers AffyCompatible).  |
| directory   | (Optional) length 1 character vector providing the full path to the local directory where the NetAffx database and files will be managed. Defaults to a session-specific temporary directory. |
| ...         | Additional arguments, not currently used.   |

**Details**

NetAffxResource creates a [NetAffxResource](#) object that manages a data base of NetAffx resources. The data base is located on the user file system as indicated by the directory argument. It is populated when the user requests specific files.

Users should provide a directory argument, so that the data base can be reused across sessions. This means that large files are only downloaded once.

Additional details are provided in the vignette accompanying this package.

**Value**

[NetAffxResource](#).

**Author(s)**

Martin Morgan

**See Also**

[NetAffxResource](#). The vignette 'Annotations with NetAffx'.

---

NetAffxResource-class *Manage NetAffx annotations*

---

### Description

This class is for accessing and retrieving NetAffx annotation files for Affymetrix chips. It requires that the user has established an account with Affymetrix. Detailed usage is described in the 'Annotations with NetAffx' vignette.

### Objects from the Class

Objects are created by calls of the form `NetAffxResource(user="user@research.org", password="terces")`, as illustrated below and in the 'Annotations with NetAffx vignette'.

### Slots

**user:** Object of class "character" containing the full user name (typically an email address) obtained when the user registers with Affymetrix.

**password:** Object of class "character" containing the password associated with the Affymetrix user name. The password is entered and stored in plain text, and should not be considered secure within Bioconductor.

**affxLicence:** Object of class "character" containing the licence, obtained from Affymetrix, for retrieving NetAffx annotations. The licence is usually software specific (e.g., for the Affy-Compatible package of Bioconductor). Users will not usually change the default value.

**directory:** Object of class "character", containing the local directory path where the NetAffx data base and annotation files will reside.

**annotationsFile:** Object of class "character" containing the file name of the NetAffx data base. Users will not usually change the default value.

**affxUrl:** Object of class "character" containing the URL used to download the NetAffx annotations file. Users will not usually change the default value.

### Extends

Class "[NetAffxCompatibility](#)", directly. Class "[AffyCompatible](#)", by class "NetAffxCompatibility", distance 2.

### Methods

[[ signature(x = "NetAffxResource", i = "numeric", j = "missing")

[[ signature(x = "NetAffxResource", i = "character", j = "missing"): Retrieve an object representing all NetAffx annotations available for an array. The argument *i* may be a numeric index or character string corresponding to an array name returned by `names(x)`.

[[ signature(x = "NetAffxResource", i = "numeric", j = "numeric")

[[ signature(x = "NetAffxResource", i = "character", j = "numeric")

[[ signature(x = "NetAffxResource", i = "numeric", j = "character")

**[[ signature(x = "NetAffxResource", i = "character", j = "character"):** Retrieve an object representing a specific annotation for an array. The argument *i* may be a numeric index or character string corresponding to an array name returned by `names(x)`. The argument *j* may be a numeric index or character string corresponding to an annotation returned by `description(x[[i]])`.

**affxLicence** signature(x = "NetAffxResource"): the character string representation of the licence used to access the NetAffx site.

**affxUrl** signature(x = "NetAffxResource"): the url used to access the NetAffx site.

**annotationsFile** signature(x = "NetAffxResource"): the file name of NetAffx data base, as stored on the local disk of the user.

**directory** signature(x = "NetAffxResource"): the directory in which the NetAffx data base and annotation files are stored on the local disk of the user.

**names** signature(x = "NetAffxResource"): the names of the arrays for which annotations are available.

**password** signature(x = "NetAffxResource"): The password associated with the user name obtained when the user registered with Affymetrix. The password is entered and stored as plain text, and is not secure.

**readAnnotation** signature(netAffxResource = "NetAffxResource", array = "numeric", annotation = "numeric")

**readAnnotation** signature(netAffxResource = "NetAffxResource", array = "numeric", annotation = "character")

**readAnnotation** signature(netAffxResource = "NetAffxResource", array = "character", annotation = "numeric")

**readAnnotation** signature(netAffxResource = "NetAffxResource", array = "character", annotation = "character")

**readAnnotation** signature(netAffxResource = "NetAffxResource", array = "missing", annotation = "AffxAnnotation")

Read a specific annotation file as an R object. For the first four methods, *array*, if present, indexes or names the array while *annotation* indexes or names the annotation. All methods also accept *content = "logical"* to indicate whether the content of the file should be returned if possible (*content=TRUE*, default) or only the path to the local file (*content=FALSE*) and *update = "logical"* indicating whether the file should be read from disk if available (*update=FALSE*, default) or retrieved from the NetAffx web site (*update=TRUE*). `readAnnotation` parses CSV, Tabular, and PSI files to data frames, and reads FASTA files to a list of description / sequence pairs (see `Biostrings::readFASTA`).

The `...` argument is passed to corresponding read functions. CSV files are parsed using `read.csv(conn, ...)`, Tabular files with `read.delim(conn, ...)`, PSI files with `read.delim(conn, header=FALSE, skip=1, sep="\t", ...)`, and FASTA with `readFASTA(conn, ...)`.

**show** signature(object = "NetAffxResource"): compactly show the NetAffxResource object.

**user** signature(x = "NetAffxResource"): The user name (typically email address) obtained when the user registers with the Affymetrix web site.

## Author(s)

Martin Morgan <mtmorgan@fhcrc.org>

## See Also

Vignette ‘Annotations with NetAffx’. Affymetrix user login site <https://www.affymetrix.com/site/login/login.affx>.

`.Rprofile` for ways to automatically set variables (such as password, and directory) for each session.

## Examples

```
## password must be obtained by the user
rsrc <- NetAffxResource(user="mtmorgan@fhcrc.org", password=password)
head(names(rsrc))
affxDescription(rsrc[["Bovine"]])
## paradigm to display all elements of annotation
sapply(affxAnnotation(rsrc[["Bovine"]]), force)
## obtain and read a single annotation
anno <- rsrc[["Porcine", "Annotations, CSV format"]]
df <- readAnnotation(rsrc, annotation=anno)
```

---

|         |  |
|---------|--|
| readArr | <i>Read ARR sample attribute files</i> |
|---------|--|

---

## Description

This functions reads the content of Affymetrix ARR sample attribute files created by GeneChip(R) Command Console. Objects are represented as R classes. Navigation and extraction of data are described in the vignette ‘Retrieving MAGE and ARR sample attributes’.

## Usage

```
readArr(fls, ...)
```

## Arguments

|                  |  |
|------------------|--|
| <code>fls</code> | A character vector containing the paths to files to be read. |
| <code>...</code> | Additional arguments, currently unused.                      |

## Value

An object or list of objects of class `ArrayFileSet`, depending on whether one or several file paths are provided.

## Author(s)

Martin Morgan



**See Also**

The vignette ‘Retrieving MAGE and ARR sample attributes’ for detailed illustration of use. [xclass](#) and [readXml](#) for advanced access methods.

**Examples**

```
dir <- system.file("extdata", "ARR", package="AffyCompatible")
fls <- list.files(dir, full=TRUE)
readArr(fls[[1]])
readArr(fls)
```

---

readMage

*Read DTT (MAGE) sample attribute files*

---

**Description**

This functions reads the content of Affymetrix DTT (MAGE) sample attribute files created by GCOS. Objects are represented as R classes. Navigation and extraction of data are described in the vignette ‘Retrieving MAGE and ARR sample attributes’.

**Usage**

```
readMage(fls, ...)
```

**Arguments**

`fls` A character vector containing the paths to files to be read.  
`...` Additional arguments, currently unused.

**Value**

An object or list of objects of class `ArrayFileSet`, depending on whether one or several file paths are provided.

**Author(s)**

Martin Morgan

**See Also**

The vignette ‘Retrieving MAGE and ARR sample attributes’ for detailed illustration of use. [xclass](#) and [readXml](#) for advanced access methods.

**Examples**

```
dir <- system.file("extdata", "DTT", package="AffyCompatible")
fls <- list.files(dir, full=TRUE)
basename(fls)
readArr(fls[[1]])
```

---

|             |   |
|-------------|---|
| readNetAffx | <i>Update or read a NetAffx data base</i> |
|-------------|---|

---

### Description

This function provides a way to read or update a NetAffx data base. The usual use for this is to force the data base description to be updated (by accessing the Affymetrix web site) or to retrieve the entire list of NetAffx resource data base. A more normal use case is to

### Usage

```
readNetAffx(netAffxResource, update = FALSE, ...)
```

### Arguments

|                 |  |
|-----------------|--|
| netAffxResource | An object of class <a href="#">NetAffxResource</a> containing the path to the local data base, and permissions for access the Affymetrix NetAffx resource. |
| update          | An object of logical(1) indicating whether the data base is to be updated (TRUE) or not.   |
| ...             | Additional arguments, not currently used.  |

### Value

An object representing the annotations found; see [NetAffxResource](#) for additional details.

### Author(s)

Martin Morgan

### Examples

```
## retrieve all annotations
## Not run:
example("NetAffxResource-class")
readNetAffx(rsrc)

## End(Not run)
```

---

|         |  |
|---------|--|
| readXml | <i>(Advanced) Read XML files of Affymetrix sample attributes or annotation</i> |
|---------|--|

---

## Description

This functions reads the content of Affymetrix DTT (MAGE) or ARR sample files, or AffxNetAffx-AnnotFile array annotation files, return an internal representation of the underlying XML. The return value can be used with [xclass](#) to rapidly create R objects representing deeply nested nodes of the document. The return value can also be used with `xpathApply` in the XML package for access to character-based representation of the data.

## Usage

```
readXml(f1, ...)
```

## Arguments

|                  |  |
|------------------|--|
| <code>f1</code>  | A length-1 character vector giving the full path to the file to be parsed. |
| <code>...</code> | Additional arguments, passed to <code>xmlTreeParse</code> .                |

## Value

An object of class `XMLInternalDocument`.

## Author(s)

Martin Morgan

## See Also

The vignettes ‘Retrieving MAGE and ARR sample attributes’ and ‘Annotation retrieval with NetAffxResource’ for detailed illustration of use.

[xclass](#) for instantiating objects from xpath queries.

## Examples

```
mageDir <- system.file("extdata", "DTT", package="AffyCompatible")
mageFiles <- list.files(mageDir, full=TRUE)
xml <- readXml(mageFiles[[1]])
xclass(xml, "//MeasuredBioAssay")[[1]]
```

---

`xclass`*(Advanced) Construct R objects from XML document queries*

---

**Description**

This function takes an xml object, created from [readXml](#), and an xpath query referencing an element of the document, and returns an R object representing the element.

The xpath query must retrieve an element, not an attribute.

**Usage**

```
xclass(doc, xpathq, prefix = character(0))
```

**Arguments**

|                     |  |
|---------------------|--|
| <code>doc</code>    | An XMLInternalDocument retrieved using <a href="#">readXml</a> .   |
| <code>xpathq</code> | A length-1 character string representing an xpath query into the document.   |
| <code>prefix</code> | A prefix used to map, when necessary, between the XML class name and the corresponding R class. For NetAffx-based classes, the prefix is 'Affx'; no prefix is required for DTT or ARR documents. |

**Value**

An R object corresponding to the class(es) referenced by the xpath query.

**Author(s)**

Martin Morgan

**References**

<http://www.w3.org/TR/xpath> provides a very useful description of xpath. Additional references are in the vignette 'Retrieving MAGE and ARR sample attributes'.

**See Also**

The vignettes 'Retrieving MAGE and ARR sample attributes' and 'Annotation retrieval with NetAffxResource' for detailed illustration of use.

[readXml](#) for appropriately reading XML documents.

**Examples**

```
example(readXml)
```

---

xmlValue.XMLAttributeValue

*Additional functions defined in AffyCompatible*


---

## Description

This page documents additional, miscellaneous, functions defined by AffyCompatible

## Usage

```
## S3 method for class XMLAttributeValue
xmlValue(x, ignoreComments = FALSE, recursive = TRUE,
         encoding=XML:::CE_NATIVE, trim=FALSE)
```

## Arguments

|                |   |
|----------------|---|
| x              | An object of class XMLAttributeValue.                     |
| ignoreComments | Logical; ignored by the xmlValue.XMLAttributeValue method |
| recursive      | ignored   |
| encoding       | File encoding   |
| trim           | Remove trailing whitespace? Ignored by this method        |

## Value

A character vector (usually length 1) containing the character representation of the XML attribute.

---

zzz-ARRCompatibility-class

*(Auto-generated) classes and methods for navigating ARR sample attributes.*


---

## Description

These classes and methods facilitate navigation of the Affymetrix ARR sample attribute files. ARR files are generated by the Affymetrix AGCC (GeneChip(R) Command Console) application. Specifically, the classes are R representations of the Affymetrix classes defined in ArraySetAndTemplate-File.dtd. The 'show' methods for these classes indicate how objects are to be accessed; details are in the 'Retrieving MAGe and ARR sample attributes' vignette.

## Objects from the Class

Objects are created internally, by parsing XML retrieved from ARR, and are not created directly by the user.

**Slots**

Objects have a variety of slots. To discover slot names and content type, see the `slotNames` portion of the example, below.

**Extends**

Class "[ARRCompatibility](#)", directly. Class "[AffyCompatible](#)", by class "ARRCompatibility", distance 2.

**Methods**

'Accessor' methods are defined on auto-generated classes to access slot content.

Accessors are apparent from displaying the object; each line of the display is of the form `accessor: content`. To retrieve the content of an instance object, invoke `accessor(object)`.

Some objects contains several instances of another object. These 'typed associations' are represented by the [.TypedAssociation](#) class, and are navigated or subset as one would a standard R vector, e.g., using `[, []`. A useful paradigm for displaying all instances is illustrated in the vignette.

**Author(s)**

Martin Morgan <[mtmorgan@fhcrc.org](mailto:mtmorgan@fhcrc.org)>

**See Also**

Vignette 'Retrieving MAGE and ARR sample attributes'.

**Examples**

```
## defined and auto-generated classes
names(slot(getClass("ARRCompatibility",
                  where=getNamespace("AffyCompatible")),
       "subclasses"))
```

---

zzz-DTTCompatibility-class

*(Auto-generated) classes and methods for navigating DTT sample attributes.*

---

**Description**

These classes and methods facilitate navigation of the Affymetrix DTT sample attribute files. DTT files are generated by the Affymetrix GCOS application. Specifically, the classes are R representations of the Affymetrix classes defined in MAGE-ML.dtd. The 'show' methods for these classes indicate how objects are to be accessed; details are in the 'Retrieving MAGE and ARR sample attributes' vignette.

## Objects from the Class

Objects are created internally, by parsing XML retrieved from ARR, and are not created directly by the user.

## Slots

Objects have a variety of slots. To discover slot names and content type, see the `slotNames` portion of the example, below.

## Extends

Class "[DTTCompatibility](#)", directly. Class "[AffyCompatible](#)", by class "DTTCompatibility", distance 2.

## Methods

'Accessor' methods are defined on auto-generated classes to access slot content.

Accessors are apparent from displaying the object; each line of the display is of the form `accessor: content`. To retrieve the content of an instance object, invoke `accessor(object)`.

Some objects contains several instances of another object. These 'typed associations' are represented by the [.TypedAssociation](#) class, and are navigated or subset as one would a standard R vector, e.g., using `[, []`. A useful paradigm for displaying all instances is illustrated in the vignette.

## Author(s)

Martin Morgan <[mtmorgan@fhcrc.org](mailto:mtmorgan@fhcrc.org)>

## See Also

Vignette 'Retrieving MAGE and ARR sample attributes'.

## Examples

```
## defined and auto-generated classes
names(slot(getClass("DTTCompatibility",
                  where=getNamespace("AffyCompatible")),
       "subclasses"))
```

---

zzz-NetAffxCompatibility-class

*(Auto-generated) classes and methods for navigating NetAffx resources*

---

## Description

These classes and methods facilitate navigation of the NetAffx resource data base. They are R representations of the Affymetrix classes defined in NetAffxAnnotFileList.dtd. The 'show' methods for these classes indicate how objects are to be accessed; details are in the 'Annotations with NetAffx' vignette.

The auto-generated classes are discoverable as indicated in the example below.

## Objects from the Class

Objects are created internally, by parsing XML retrieved from the local NetAffx data base or from the Affymetrix web site.

## Slots

Objects have a variety of slots. To discover slot names and content type, see the slotNames portion of the example, below.

## Extends

Class "[NetAffxCompatibility](#)", directly. Class "[AffyCompatible](#)", by class "[NetAffxCompatibility](#)", distance 2.

## Methods

'Read' methods are defined on auto-generated classes to access slot content.

Accessors are apparent from displaying the object; each line of the display is of the form accessor: content. To retrieve the content of an instance object, invoke accessor(object).

Some objects contains several instances of another object. These 'typed associations' are represented by the [.TypedAssociation](#) class, and are navigated or subset as one would a standard R vector, e.g., using [, [[]]. A useful paradigm for displaying all instances is illustrated on the [NetAffxResource](#) page.

## Author(s)

Martin Morgan <mtmorgan@fhcrc.org>



### See Also

Vignette 'Annotations with NetAffx'. Affymetrix user login site <https://www.affymetrix.com/site/login/login.affx>.

[NetAffxResource](#) for creating and managing the NetAffx data base.

[NetAffxResource](#) for navigating the data base.

### Examples

```
## defined and auto-generated classes
getClass("NetAffxCompatibility", where=getNamespace("AffyCompatible"))

## slots of an auto-generated class
getSlots(getClass("AffxNetAffxAnnotFileList",
  where=getNamespace("AffyCompatible")))

## Not run:
example("NetAffxrResource-class")

## End(Not run)
```

# Index

- \*Topic **classes**
  - AffyCompatible-class, 3
  - Internal-class, 4
  - NetAffxResource-class, 6
  - readArr, 8
  - readMage, 9
  - zzz-ARRCompatibility-class, 13
  - zzz-DTTCompatibility-class, 14
  - zzz-NetAffxCompatibility-class, 16
- \*Topic **database**
  - NetAffxResource, 4
  - readNetAffx, 10
- \*Topic **manip**
  - readXml, 11
  - xclass, 12
  - xmlValue.XMLAttributeValue, 13
- \*Topic **methods**
  - accessors, 2
- \*Topic **package**
  - AffyCompatible-package, 2
  - .Rprofile, 8
  - .TypedAssociation, 14–16
  - .TypedAssociation-class
    - (Internal-class), 4
  - [, .TypedAssociation, ANY, ANY-method
    - (Internal-class), 4
  - [, .TypedAssociation, numeric, missing, ANY-method
    - (Internal-class), 4
  - [, .TypedAssociation, numeric, missing-method
    - (Internal-class), 4
  - [[, NetAffxResource, character, character-method
    - (NetAffxResource-class), 6
  - [[, NetAffxResource, character, missing-method
    - (NetAffxResource-class), 6
  - [[, NetAffxResource, character, numeric-method
    - (NetAffxResource-class), 6
  - [[, NetAffxResource, numeric, character-method
    - (NetAffxResource-class), 6
  - [[, NetAffxResource, numeric, missing-method
    - (NetAffxResource-class), 6
  - [[, NetAffxResource, numeric, numeric-method
    - (NetAffxResource-class), 6
  - accession (zzz-DTTCompatibility-class), 14
  - accession, DatabaseEntry-method
    - (zzz-DTTCompatibility-class), 14
  - accessions\_assnlist
    - (zzz-DTTCompatibility-class), 14
  - accessions\_assnlist, BibliographicReference-method
    - (zzz-DTTCompatibility-class), 14
  - Accessions\_assnlist-class
    - (zzz-DTTCompatibility-class), 14
  - accessionVersion
    - (zzz-DTTCompatibility-class), 14
  - accessionVersion, DatabaseEntry-method
    - (zzz-DTTCompatibility-class), 14
  - accessors, 2
  - acpassword (NetAffxResource), 4
  - action (zzz-DTTCompatibility-class), 14
  - action, Audit-method
    - (zzz-DTTCompatibility-class), 14
  - action\_assn
    - (zzz-DTTCompatibility-class), 14
  - action\_assn, Treatment-method
    - (zzz-DTTCompatibility-class), 14
  - Action\_assn-class
    - (zzz-DTTCompatibility-class), 14
  - actionMeasurement\_assn

- (zzz-DTTCompatibility-class),  
14
- actionMeasurement\_assn,Treatment-method  
(zzz-DTTCompatibility-class),  
14
- ActionMeasurement\_assn-class  
(zzz-DTTCompatibility-class),  
14
- activityDate  
(zzz-DTTCompatibility-class),  
14
- activityDate,ProtocolApplication-method  
(zzz-DTTCompatibility-class),  
14
- address(zzz-DTTCompatibility-class), 14
- address,Organization-method  
(zzz-DTTCompatibility-class),  
14
- address,Person-method  
(zzz-DTTCompatibility-class),  
14
- adjustments\_assnlist  
(zzz-DTTCompatibility-class),  
14
- adjustments\_assnlist,ArrayManufactureDeviationFileRecord  
(zzz-DTTCompatibility-class),  
14
- Adjustments\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- affiliation\_assnref  
(zzz-DTTCompatibility-class),  
14
- affiliation\_assnref,Person-method  
(zzz-DTTCompatibility-class),  
14
- Affiliation\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- affxAnnotation  
(zzz-NetAffxCompatibility-class),  
16
- affxAnnotation,AffxArray-method  
(zzz-NetAffxCompatibility-class),  
16
- AffxAnnotation-class  
(zzz-NetAffxCompatibility-class),  
16
- affxArray  
(zzz-NetAffxCompatibility-class),  
16
- affxArray,AffxNetAffxAnnotFileList-method  
(zzz-NetAffxCompatibility-class),  
16
- affxArray,Array\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- AffxArray-class  
(zzz-NetAffxCompatibility-class),  
16
- affxCatgory  
(zzz-DTTCompatibility-class),  
14
- affxCatgory,OntologyEntry-method  
(zzz-DTTCompatibility-class),  
14
- affxComment  
(zzz-NetAffxCompatibility-class),  
16
- affxComment,AffxAnnotation-method  
(zzz-NetAffxCompatibility-class),  
16
- affxComment,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- AffxComment-class  
(zzz-NetAffxCompatibility-class),  
16
- affxCompression  
(zzz-NetAffxCompatibility-class),  
16
- affxCompression,AffxUrl-method  
(zzz-NetAffxCompatibility-class),  
16
- affxCrc32  
(zzz-NetAffxCompatibility-class),  
16
- affxCrc32,AffxFile-method  
(zzz-NetAffxCompatibility-class),  
16
- affxCrc32,AffxUrl-method  
(zzz-NetAffxCompatibility-class),  
16
- affxDate  
(zzz-NetAffxCompatibility-class),  
16

- affxDate, AffxFile-method  
 (zzz-NetAffxCompatibility-class),  
 16
- affxDate, Audit-method  
 (zzz-DTTCompatibility-class),  
 14
- affxDescription  
 (zzz-NetAffxCompatibility-class),  
 16
- affxDescription, AffxAnnotation-method  
 (zzz-NetAffxCompatibility-class),  
 16
- affxDescription, AffxArray-method  
 (zzz-NetAffxCompatibility-class),  
 16
- affxEnd (zzz-DTTCompatibility-class), 14
- affxEnd, CompositePosition-method  
 (zzz-DTTCompatibility-class),  
 14
- affxEnd, ReporterPosition-method  
 (zzz-DTTCompatibility-class),  
 14
- affxEnd, SequencePosition-method  
 (zzz-DTTCompatibility-class),  
 14
- affxFile  
 (zzz-NetAffxCompatibility-class),  
 16
- affxFile, AffxAnnotation-method  
 (zzz-NetAffxCompatibility-class),  
 16
- AffxFile-class  
 (zzz-NetAffxCompatibility-class),  
 16
- affxImage (zzz-DTTCompatibility-class),  
 14
- affxImage, PhysicalBioAssayData\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- affxLicence  
 (zzz-NetAffxCompatibility-class),  
 16
- affxLicence, NetAffxResource-method  
 (NetAffxResource-class), 6
- affxName  
 (zzz-NetAffxCompatibility-class),  
 16
- affxName, AffxArray-method  
 (zzz-NetAffxCompatibility-class),  
 16
- affxName, AffxFile-method  
 (zzz-NetAffxCompatibility-class),  
 16
- AffxNetAffxAnnotFileList-class  
 (zzz-NetAffxCompatibility-class),  
 16
- affxOrder (zzz-DTTCompatibility-class),  
 14
- affxOrder, BioDataCube-method  
 (zzz-DTTCompatibility-class),  
 14
- affxOrder, Treatment-method  
 (zzz-DTTCompatibility-class),  
 14
- affxPerson  
 (zzz-DTTCompatibility-class),  
 14
- affxPerson, Contact\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- affxPerson, Person\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- affxRow (zzz-DTTCompatibility-class), 14
- affxRow, FeatureLocation-method  
 (zzz-DTTCompatibility-class),  
 14
- affxRow, Zone-method  
 (zzz-DTTCompatibility-class),  
 14
- affxSequence  
 (zzz-DTTCompatibility-class),  
 14
- affxSequence, BioSequence-method  
 (zzz-DTTCompatibility-class),  
 14
- affxSize  
 (zzz-NetAffxCompatibility-class),  
 16
- affxSize, AffxFile-method  
 (zzz-NetAffxCompatibility-class),  
 16
- affxSize, AffxUrl-method  
 (zzz-NetAffxCompatibility-class),  
 16
- affxStart (zzz-DTTCompatibility-class),

- 14
- affxStart, CompositePosition-method  
(zzz-DTTCompatibility-class),  
14
- affxStart, ReporterPosition-method  
(zzz-DTTCompatibility-class),  
14
- affxStart, SequencePosition-method  
(zzz-DTTCompatibility-class),  
14
- affxStatus  
(zzz-NetAffxCompatibility-class),  
16
- affxStatus, AffxNetAffxAnnotFileList-method  
(zzz-NetAffxCompatibility-class),  
16
- AffxStatus-class  
(zzz-NetAffxCompatibility-class),  
16
- affxText (zzz-DTTCompatibility-class),  
14
- affxText, Description-method  
(zzz-DTTCompatibility-class),  
14
- affxText, Protocol-method  
(zzz-DTTCompatibility-class),  
14
- affxTitle (zzz-DTTCompatibility-class),  
14
- affxTitle, BibliographicReference-method  
(zzz-DTTCompatibility-class),  
14
- affxTitle, Protocol-method  
(zzz-DTTCompatibility-class),  
14
- affxType  
(zzz-NetAffxCompatibility-class),  
16
- affxType, AffxAnnotation-method  
(zzz-NetAffxCompatibility-class),  
16
- affxType, AffxFile-method  
(zzz-NetAffxCompatibility-class),  
16
- affxUrl  
(zzz-NetAffxCompatibility-class),  
16
- affxUrl, AffxFile-method  
(zzz-NetAffxCompatibility-class),  
16
- affxUrl, NetAffxResource-method  
(NetAffxResource-class), 6
- AffxUrl-class  
(zzz-NetAffxCompatibility-class),  
16
- affyBarcode  
(zzz-ARRCompatibility-class),  
13
- affyBarcode, PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- AffyCompatible, 6, 14–16
- AffyCompatible  
(AffyCompatible-package), 2
- AffyCompatible-class, 3
- AffyCompatible-package, 2
- analysisResults\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- analysisResults\_assnreflist, Experiment-method  
(zzz-DTTCompatibility-class),  
14
- AnalysisResults\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- annotations\_assnlist  
(zzz-DTTCompatibility-class),  
14
- annotations\_assnlist, Description-method  
(zzz-DTTCompatibility-class),  
14
- annotations\_assnlist, ExperimentalFactor-method  
(zzz-DTTCompatibility-class),  
14
- Annotations\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- annotationsFile (accessors), 2
- annotationsFile, NetAffxResource-method  
(NetAffxResource-class), 6
- Array-class  
(zzz-DTTCompatibility-class),  
14
- array\_assnlist  
(zzz-DTTCompatibility-class),  
14

- array\_assnlist,Array\_package-method  
(zzz-DTTCompatibility-class),  
[14](#)
- Array\_assnlist-class  
(zzz-DTTCompatibility-class),  
[14](#)
- array\_assnref  
(zzz-DTTCompatibility-class),  
[14](#)
- array\_assnref,BioAssayCreation-method  
(zzz-DTTCompatibility-class),  
[14](#)
- array\_assnref,Hybridization-method  
(zzz-DTTCompatibility-class),  
[14](#)
- Array\_assnref-class  
(zzz-DTTCompatibility-class),  
[14](#)
- array\_package  
(zzz-DTTCompatibility-class),  
[14](#)
- array\_package,MAGE-ML-method  
(zzz-DTTCompatibility-class),  
[14](#)
- Array\_package-class  
(zzz-DTTCompatibility-class),  
[14](#)
- array\_ref (zzz-DTTCompatibility-class),  
[14](#)
- array\_ref,Array\_assnref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- array\_ref,Arrays\_assnreflist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- Array\_ref-class  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayAttribute  
(zzz-ARRCompatibility-class),  
[13](#)
- arrayAttribute,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
[13](#)
- ArrayAttribute-class  
(zzz-ARRCompatibility-class),  
[13](#)
- arrayDesign  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayDesign,ArrayDesign\_assnlist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ArrayDesign-class  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayDesign\_assnlist  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayDesign\_assnlist,ArrayDesign\_package-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ArrayDesign\_assnlist-class  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayDesign\_assnref  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayDesign\_assnref,Array-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ArrayDesign\_assnref-class  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayDesign\_package  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayDesign\_package,MAGE-ML-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ArrayDesign\_package-class  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayDesign\_ref  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayDesign\_ref,ArrayDesign\_assnref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ArrayDesign\_ref-class  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayGroup  
(zzz-DTTCompatibility-class),  
[14](#)
- arrayGroup,ArrayGroup\_assnlist-method

- (zzz-DTTCompatibility-class),  
14
- ArrayGroup-class  
(zzz-DTTCompatibility-class),  
14
- arrayGroup\_assnlist  
(zzz-DTTCompatibility-class),  
14
- arrayGroup\_assnlist,Array\_package-method  
(zzz-DTTCompatibility-class),  
14
- ArrayGroup\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- arrayGroup\_assnref  
(zzz-DTTCompatibility-class),  
14
- arrayGroup\_assnref,Array-method  
(zzz-DTTCompatibility-class),  
14
- ArrayGroup\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- arrayGroup\_ref  
(zzz-DTTCompatibility-class),  
14
- arrayGroup\_ref,ArrayGroup\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- ArrayGroup\_ref-class  
(zzz-DTTCompatibility-class),  
14
- arrayIdentifier  
(zzz-DTTCompatibility-class),  
14
- arrayIdentifier,Array-method  
(zzz-DTTCompatibility-class),  
14
- arrayManufacture  
(zzz-DTTCompatibility-class),  
14
- arrayManufacture,ArrayManufacture\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ArrayManufacture-class  
(zzz-DTTCompatibility-class),  
14
- arrayManufacture\_assnlist  
(zzz-DTTCompatibility-class),  
14
- arrayManufacture\_assnlist,Array\_package-method  
(zzz-DTTCompatibility-class),  
14
- ArrayManufacture\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- arrayManufacture\_ref  
(zzz-DTTCompatibility-class),  
14
- arrayManufacture\_ref,Information\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- ArrayManufacture\_ref-class  
(zzz-DTTCompatibility-class),  
14
- arrayManufactureDeviation  
(zzz-DTTCompatibility-class),  
14
- arrayManufactureDeviation,ArrayManufactureDeviations\_assnlist  
(zzz-DTTCompatibility-class),  
14
- ArrayManufactureDeviation-class  
(zzz-DTTCompatibility-class),  
14
- arrayManufactureDeviations\_assnlist  
(zzz-DTTCompatibility-class),  
14
- arrayManufactureDeviations\_assnlist,Array-method  
(zzz-DTTCompatibility-class),  
14
- ArrayManufactureDeviations\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- arrayManufacturers\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- arrayManufacturers\_assnreflist,ArrayManufacture-method  
(zzz-DTTCompatibility-class),  
14
- ArrayManufacturers\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- arrayName (zzz-ARRCompatibility-class),  
13
- arrayName,PhysicalArray-method  
(zzz-ARRCompatibility-class),

- 13
- arrays\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- arrays\_assnreflist,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- arrays\_assnreflist,ArrayManufacture-method  
(zzz-DTTCompatibility-class),  
14
- Arrays\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- ArraySetFile-class  
(zzz-ARRCompatibility-class),  
13
- arraySpacingX  
(zzz-DTTCompatibility-class),  
14
- arraySpacingX,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- arraySpacingY  
(zzz-DTTCompatibility-class),  
14
- arraySpacingY,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- arrayXOrigin  
(zzz-DTTCompatibility-class),  
14
- arrayXOrigin,Array-method  
(zzz-DTTCompatibility-class),  
14
- arrayYOrigin  
(zzz-DTTCompatibility-class),  
14
- arrayYOrigin,Array-method  
(zzz-DTTCompatibility-class),  
14
- ARRCompatibility, 14
- ARRCompatibility-class  
(AffyCompatible-class), 3
- associations\_assnlist  
(zzz-DTTCompatibility-class),  
14
- associations\_assnlist,OntologyEntry-method  
(zzz-DTTCompatibility-class),  
14
- Associations\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- audit (zzz-DTTCompatibility-class), 14
- audit,AuditTrail\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Audit-class  
(zzz-DTTCompatibility-class),  
14
- auditAndSecurity\_package  
(zzz-DTTCompatibility-class),  
14
- auditAndSecurity\_package,MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- AuditAndSecurity\_package-class  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Array-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ArrayManufacture-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Audit-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,BibliographicReference-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,BioAssayCreation-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,BioAssayDataCluster-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,BioAssayDimension-method



- (zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,BioAssayMap-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,BioAssayTreatment-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,BioSample-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,BioSequence-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,BioSource-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Channel-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,CompositeCompositeMap-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,CompositeGroup-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,CompositeSequence-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,CompositeSequenceDimensionalMethod  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Compound-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Database-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,DerivedBioAssay-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,DerivedBioAssayData-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,DerivedSignal-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Description-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Error-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ExpectedValue-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Experiment-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ExperimentalFactor-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ExperimentDesign-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,FactorValue-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Failed-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Feature-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,FeatureDimension-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,FeatureExtraction-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,FeatureGroup-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,FeatureReporterMap-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Fiducial-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Hardware-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,HardwareApplication-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Hybridization-method  
(zzz-DTTCompatibility-class),  
14

- (zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Image-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ImageAcquisition-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,LabeledExtract-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ManufactureLIMS-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ManufactureLIMSBiomaterials-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,MeasuredBioAssay-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,MeasuredBioAssayData-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,MeasuredSignal-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Node-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,NodeContents-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Organization-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Parameter-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Person-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,PhysicalArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,PhysicalBioAssay-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,PresentAbsent-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Protocol-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ProtocolApplication-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,PValue-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,QuantitationTypeDimension-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,QuantitationTypeMap-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Ratio-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Reporter-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ReporterCompositeMap-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ReporterDimension-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,ReporterGroup-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Security-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,SecurityGroup-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,SeqFeature-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,Software-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist,SoftwareApplication-method  
(zzz-DTTCompatibility-class),  
14

- (zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist, SpecializedQuantitationTypeAssayAssnlist-class, BioAssay\_package-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist, Transformation-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist, Treatment-method  
(zzz-DTTCompatibility-class),  
14
- auditTrail\_assnlist, Zone-method  
(zzz-DTTCompatibility-class),  
14
- AuditTrail\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- authors (zzz-DTTCompatibility-class), 14
- authors, BibliographicReference-method  
(zzz-DTTCompatibility-class),  
14
- barcode (zzz-DTTCompatibility-class), 14
- barcode, ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- basis (zzz-DTTCompatibility-class), 14
- basis, SeqFeature-method  
(zzz-DTTCompatibility-class),  
14
- bibliographicReference  
(zzz-DTTCompatibility-class),  
14
- bibliographicReference, BibliographicReferences\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- BibliographicReference-class  
(zzz-DTTCompatibility-class),  
14
- bibliographicReferences\_assnlist  
(zzz-DTTCompatibility-class),  
14
- bibliographicReferences\_assnlist, DescriptionBioAssayCreation-assn-method  
(zzz-DTTCompatibility-class),  
14
- BibliographicReferences\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssay\_assnlist  
(zzz-DTTCompatibility-class),  
14
- bioAssay\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssay\_assnref  
(zzz-DTTCompatibility-class),  
14
- bioAssay\_assnref, BioAssayTuple-method  
(zzz-DTTCompatibility-class),  
14
- BioAssay\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- bioAssay\_package  
(zzz-DTTCompatibility-class),  
14
- bioAssay\_package, MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- BioAssay\_package-class  
(zzz-DTTCompatibility-class),  
14
- BioAssay\_ref-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayCreation  
(zzz-DTTCompatibility-class),  
14
- bioAssayCreation, BioAssayCreation\_assn-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayCreation-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayCreation\_assn  
(zzz-DTTCompatibility-class),  
14
- bioAssayCreation\_assn, PhysicalBioAssay-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayCreation\_assn-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayData\_assnlist

- (zzz-DTTCompatibility-class),  
14
- bioAssayData\_assnlist, BioAssayData\_package-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayData\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayData\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- bioAssayData\_assnreflist, Experiment-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayData\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayData\_package  
(zzz-DTTCompatibility-class),  
14
- bioAssayData\_package, MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayData\_package-class  
(zzz-DTTCompatibility-class),  
14
- BioAssayData\_ref-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayDataCluster  
(zzz-DTTCompatibility-class),  
14
- bioAssayDataCluster, BioAssayDataCluster\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayDataCluster-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayDataCluster\_assnlist  
(zzz-DTTCompatibility-class),  
14
- bioAssayDataCluster\_assnlist, HigherLevelAnalysis-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayDataCluster\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayDataCluster\_ref  
(zzz-DTTCompatibility-class),  
14
- bioAssayDataCluster\_ref, AnalysisResults\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayDataCluster\_ref-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayDataSources\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- bioAssayDataSources\_assnreflist, Transformation-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayDataSources\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayDimension  
(zzz-DTTCompatibility-class),  
14
- bioAssayDimension, BioAssayDimension\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayDimension-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayDimension\_assnlist  
(zzz-DTTCompatibility-class),  
14
- bioAssayDimension\_assnlist, BioAssayData\_package-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayDimension\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayDimension\_assnref  
(zzz-DTTCompatibility-class),  
14
- bioAssayDimension\_assnref, DerivedBioAssayData-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayDimension\_assnref, MeasuredBioAssayData-method  
(zzz-DTTCompatibility-class),  
14
- bioAssayDimension\_assnref, NodeContents-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayDimension\_assnref-class

- (zzz-DTTCompatibility-class),  
14
- bioAssayDimension\_ref  
(zzz-DTTCompatibility-class),  
14
- bioAssayDimension\_ref,BioAssayDimension\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayDimension\_ref-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayFactorValues\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- bioAssayFactorValues\_assnreflist,DerivedBioAssayFactorValues-assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- bioAssayFactorValues\_assnreflist,MeasuredBioAssayFactorValues-assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- bioAssayFactorValues\_assnreflist,PhysicalBioAssayFactorValues-assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayFactorValues\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayMap  
(zzz-DTTCompatibility-class),  
14
- bioAssayMap,BioAssayMap\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayMap-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayMap\_assnlist  
(zzz-DTTCompatibility-class),  
14
- bioAssayMap\_assnlist,BioAssayData\_package-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayMap\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayMap\_ref  
(zzz-DTTCompatibility-class),  
14
- bioAssayMap\_ref,BioAssayMaps\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- (zzz-DTTCompatibility-class),  
14
- bioAssayMap\_ref,DerivedBioAssayMap\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayMap\_ref-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayMapping  
(zzz-DTTCompatibility-class),  
14
- bioAssayMapping,BioAssayMapping\_assn-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayMapping-class  
(zzz-DTTCompatibility-class),  
14
- BioAssayMapping\_assn  
(zzz-DTTCompatibility-class),  
14
- BioAssayMapping\_assn,Transformation-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayMapping\_assn-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayMaps\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- bioAssayMaps\_assnreflist,BioAssayMapping-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayMaps\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayMapTarget\_assnref  
(zzz-DTTCompatibility-class),  
14
- bioAssayMapTarget\_assnref,BioAssayMap-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayMapTarget\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- bioAssays\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- bioAssays\_assnreflist,BioAssayDimension-method  
(zzz-DTTCompatibility-class),  
14

- (zzz-DTTCompatibility-class),  
14
- bioAssays\_assnreflist, Experiment-method  
(zzz-DTTCompatibility-class),  
14
- BioAssays\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayTreatment  
(zzz-DTTCompatibility-class),  
14
- bioAssayTreatment, BioAssayTreatments\_assnlist  
(zzz-DTTCompatibility-class),  
14
- BioAssayTreatment-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayTreatments\_assnlist  
(zzz-DTTCompatibility-class),  
14
- bioAssayTreatments\_assnlist, PhysicalBioAssayTreatment  
(zzz-DTTCompatibility-class),  
14
- BioAssayTreatments\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayTuple  
(zzz-DTTCompatibility-class),  
14
- bioAssayTuple, BioAssayTuples\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayTuple-class  
(zzz-DTTCompatibility-class),  
14
- bioAssayTuples\_assnlist  
(zzz-DTTCompatibility-class),  
14
- bioAssayTuples\_assnlist, BioDataTuples-method  
(zzz-DTTCompatibility-class),  
14
- BioAssayTuples\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- bioDataCube  
(zzz-DTTCompatibility-class),  
14
- bioDataCube, BioDataValues\_assn-method  
(zzz-DTTCompatibility-class),  
14
- BioDataCube-class  
(zzz-DTTCompatibility-class),  
14
- bioDataTuples  
(zzz-DTTCompatibility-class),  
14
- bioDataTuples, BioDataValues\_assn-method  
(zzz-DTTCompatibility-class),  
14
- BioDataTuples-class  
(zzz-DTTCompatibility-class),  
14
- bioDataValues\_assn  
(zzz-DTTCompatibility-class),  
14
- bioDataValues\_assn, DerivedBioAssayData-method  
(zzz-DTTCompatibility-class),  
14
- bioDataValues\_assn, MeasuredBioAssayData-method  
(zzz-DTTCompatibility-class),  
14
- BioDataValues\_assn-class  
(zzz-DTTCompatibility-class),  
14
- bioEvent\_package  
(zzz-DTTCompatibility-class),  
14
- bioEvent\_package, MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- BioEvent\_package-class  
(zzz-DTTCompatibility-class),  
14
- biologicalCharacteristics\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- biologicalCharacteristics\_assnreflist, CompositeSequence-method  
(zzz-DTTCompatibility-class),  
14
- BiologicalCharacteristics\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- bioMaterial\_assnlist  
(zzz-DTTCompatibility-class),  
14
- bioMaterial\_assnlist, BioMaterial\_package-method

- (zzz-DTTCompatibility-class),  
14
- BioMaterial\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- bioMaterial\_assnref  
(zzz-DTTCompatibility-class),  
14
- bioMaterial\_assnref,BioMaterialMeasurement-method  
(zzz-DTTCompatibility-class),  
14
- bioMaterial\_assnref,ManufactureLIMS-method  
(zzz-DTTCompatibility-class),  
14
- bioMaterial\_assnref,ManufactureLIMSBiomaterial-method  
(zzz-DTTCompatibility-class),  
14
- BioMaterial\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- bioMaterial\_package  
(zzz-DTTCompatibility-class),  
14
- bioMaterial\_package,MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- BioMaterial\_package-class  
(zzz-DTTCompatibility-class),  
14
- BioMaterial\_ref-class  
(zzz-DTTCompatibility-class),  
14
- bioMaterialMeasurement  
(zzz-DTTCompatibility-class),  
14
- bioMaterialMeasurement,SourceBioMaterialMeasurements-assnlist-method  
(zzz-DTTCompatibility-class),  
14
- BioMaterialMeasurement-class  
(zzz-DTTCompatibility-class),  
14
- bioMaterialPlateCol  
(zzz-DTTCompatibility-class),  
14
- bioMaterialPlateCol,ManufactureLIMSBiomaterial-method  
(zzz-DTTCompatibility-class),  
14
- bioMaterialPlateIdentifier  
(zzz-DTTCompatibility-class),  
14
- bioMaterialPlateIdentifier,ManufactureLIMSBiomaterial-method  
(zzz-DTTCompatibility-class),  
14
- bioMaterialPlateRow  
(zzz-DTTCompatibility-class),  
14
- bioMaterialPlateRow,ManufactureLIMSBiomaterial-method  
(zzz-DTTCompatibility-class),  
14
- bioSample  
(zzz-DTTCompatibility-class),  
14
- bioSample,BioMaterial\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- BioSample-class  
(zzz-DTTCompatibility-class),  
14
- bioSample\_ref  
(zzz-DTTCompatibility-class),  
14
- bioSample\_ref,BioMaterial\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- BioSample\_ref-class  
(zzz-DTTCompatibility-class),  
14
- bioSequence  
(zzz-DTTCompatibility-class),  
14
- bioSequence,BioSequence\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- BioSequence-class  
(zzz-DTTCompatibility-class),  
14
- bioSequence\_assnlist  
(zzz-DTTCompatibility-class),  
14
- bioSequence\_assnlist,BioSequence\_package-method  
(zzz-DTTCompatibility-class),  
14
- BioSequence\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- bioSequence\_package  
(zzz-DTTCompatibility-class),

- 14
- bioSequence\_package, MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- BioSequence\_package-class  
(zzz-DTTCompatibility-class),  
14
- bioSequence\_ref  
(zzz-DTTCompatibility-class),  
14
- bioSequence\_ref, BiologicalCharacteristics\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- bioSequence\_ref, ImmobilizedCharacteristics\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- BioSequence\_ref-class  
(zzz-DTTCompatibility-class),  
14
- bioSource (zzz-DTTCompatibility-class),  
14
- bioSource, BioMaterial\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- BioSource-class  
(zzz-DTTCompatibility-class),  
14
- bioSource\_ref  
(zzz-DTTCompatibility-class),  
14
- bioSource\_ref, BioMaterial\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- BioSource\_ref-class  
(zzz-DTTCompatibility-class),  
14
- bqs\_package  
(zzz-DTTCompatibility-class),  
14
- bqs\_package, MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- BQS\_package-class  
(zzz-DTTCompatibility-class),  
14
- category\_assn  
(zzz-DTTCompatibility-class),  
14
- category\_assn, ExperimentalFactor-method  
(zzz-DTTCompatibility-class),  
14
- Category\_assn-class  
(zzz-DTTCompatibility-class),  
14
- channel (zzz-DTTCompatibility-class), 14
- channel, Channel\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- channel, Channel\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- channel, Channel\_assnlist  
(zzz-DTTCompatibility-class),  
14
- channel\_assnlist, BioAssay\_package-method  
(zzz-DTTCompatibility-class),  
14
- Channel\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- channel\_assnref  
(zzz-DTTCompatibility-class),  
14
- channel\_assnref, DerivedSignal-method  
(zzz-DTTCompatibility-class),  
14
- channel\_assnref, Error-method  
(zzz-DTTCompatibility-class),  
14
- channel\_assnref, ExpectedValue-method  
(zzz-DTTCompatibility-class),  
14
- channel\_assnref, Failed-method  
(zzz-DTTCompatibility-class),  
14
- channel\_assnref, MeasuredSignal-method  
(zzz-DTTCompatibility-class),  
14
- channel\_assnref, PresentAbsent-method  
(zzz-DTTCompatibility-class),  
14
- channel\_assnref, PValue-method  
(zzz-DTTCompatibility-class),  
14
- channel\_assnref, Ratio-method  
(zzz-DTTCompatibility-class),



- 14  
channel\_assnref, SpecializedQuantitationType-method  
(zzz-DTTCompatibility-class),  
14  
Channel\_assnref-class  
(zzz-DTTCompatibility-class),  
14  
channel\_ref  
(zzz-DTTCompatibility-class),  
14  
channel\_ref, Channel\_assnref-method  
(zzz-DTTCompatibility-class),  
14  
channel\_ref, Channels\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14  
Channel\_ref-class  
(zzz-DTTCompatibility-class),  
14  
channels\_assnreflist  
(zzz-DTTCompatibility-class),  
14  
channels\_assnreflist, DerivedBioAssay-method  
(zzz-DTTCompatibility-class),  
14  
channels\_assnreflist, Image-method  
(zzz-DTTCompatibility-class),  
14  
channels\_assnreflist, MeasuredBioAssay-method  
(zzz-DTTCompatibility-class),  
14  
channels\_assnreflist, PhysicalBioAssay-method  
(zzz-DTTCompatibility-class),  
14  
Channels\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14  
characteristics\_assnlist  
(zzz-DTTCompatibility-class),  
14  
characteristics\_assnlist, BioSample-method  
(zzz-DTTCompatibility-class),  
14  
characteristics\_assnlist, BioSource-method  
(zzz-DTTCompatibility-class),  
14  
characteristics\_assnlist, LabeledExtract-method  
(zzz-DTTCompatibility-class),  
14  
characteristics\_assnlist-class  
(zzz-DTTCompatibility-class),  
14  
clusterBioAssayData\_assnref  
(zzz-DTTCompatibility-class),  
14  
clusterBioAssayData\_assnref, BioAssayDataCluster-method  
(zzz-DTTCompatibility-class),  
14  
ClusterBioAssayData\_assnref-class  
(zzz-DTTCompatibility-class),  
14  
column (zzz-DTTCompatibility-class), 14  
column, FeatureLocation-method  
(zzz-DTTCompatibility-class),  
14  
column, Zone-method  
(zzz-DTTCompatibility-class),  
14  
componentCompounds\_assnlist  
(zzz-DTTCompatibility-class),  
14  
componentCompounds\_assnlist, Compound-method  
(zzz-DTTCompatibility-class),  
14  
ComponentCompounds\_assnlist-class  
(zzz-DTTCompatibility-class),  
14  
composite\_assnref  
(zzz-DTTCompatibility-class),  
14  
composite\_assnref, CompositePosition-method  
(zzz-DTTCompatibility-class),  
14  
Composite\_assnref-class  
(zzz-DTTCompatibility-class),  
14  
compositeCompositeMap  
(zzz-DTTCompatibility-class),  
14  
compositeCompositeMap, CompositeCompositeMap\_assnlist-method  
(zzz-DTTCompatibility-class),  
14  
compositeCompositeMap, DesignElementMap\_assnlist-method  
(zzz-DTTCompatibility-class),  
14  
CompositeCompositeMap-class

|   |   |
|---|---|
| (zzz-DTTCompatibility-class),<br>14   | (zzz-DTTCompatibility-class),<br>14   |
| compositeCompositeMap_assnlist<br>(zzz-DTTCompatibility-class),<br>14   | compositeGroup_ref<br>(zzz-DTTCompatibility-class),<br>14   |
| compositeCompositeMap_assnlist,DesignElement_package_ref,CompositeGroups_assnreflist-method<br>(zzz-DTTCompatibility-class),<br>14            | compositeGroup_ref,CompositeGroups_assnreflist-method<br>(zzz-DTTCompatibility-class),<br>14      |
| CompositeCompositeMap_assnlist-class<br>(zzz-DTTCompatibility-class),<br>14   | CompositeGroup_ref-class<br>(zzz-DTTCompatibility-class),<br>14                                   |
| compositeCompositeMap_ref<br>(zzz-DTTCompatibility-class),<br>14  | compositeGroups_assnreflist<br>(zzz-DTTCompatibility-class),<br>14                                |
| compositeCompositeMap_ref,CompositeCompositeMap_ref_assnreflist,ArrayDesign-method<br>(zzz-DTTCompatibility-class),<br>14                     | compositeGroups_assnreflist,ArrayDesign-method<br>(zzz-DTTCompatibility-class),<br>14             |
| compositeCompositeMap_ref,DesignElementMaps_assnreflist,PhysicalArrayDesign-method<br>(zzz-DTTCompatibility-class),<br>14                     | compositeGroups_assnreflist,PhysicalArrayDesign-method<br>(zzz-DTTCompatibility-class),<br>14     |
| CompositeCompositeMap_ref-class<br>(zzz-DTTCompatibility-class),<br>14  | CompositeGroups_assnreflist-class<br>(zzz-DTTCompatibility-class),<br>14                          |
| compositeCompositeMaps_assnreflist<br>(zzz-DTTCompatibility-class),<br>14   | compositePosition<br>(zzz-DTTCompatibility-class),<br>14  |
| compositeCompositeMaps_assnreflist,CompositeSequenceRef_class,CompositePositionSources_assnlist-method<br>(zzz-DTTCompatibility-class),<br>14 | compositePosition,CompositePositionSources_assnlist-method<br>(zzz-DTTCompatibility-class),<br>14 |
| CompositeCompositeMaps_assnreflist-class<br>(zzz-DTTCompatibility-class),<br>14   | compositePosition,Coordinate_assn-method<br>(zzz-DTTCompatibility-class),<br>14                   |
| compositeGroup<br>(zzz-DTTCompatibility-class),<br>14   | CompositePosition-class<br>(zzz-DTTCompatibility-class),<br>14                                    |
| compositeGroup,CompositeGroup_assnlist-method<br>(zzz-DTTCompatibility-class),<br>14  | compositePositionSources_assnlist<br>(zzz-DTTCompatibility-class),<br>14                          |
| CompositeGroup-class<br>(zzz-DTTCompatibility-class),<br>14   | compositePositionSources_assnlist,CompositeCompositeMap-me<br>(zzz-DTTCompatibility-class),<br>14 |
| compositeGroup_assnlist<br>(zzz-DTTCompatibility-class),<br>14  | CompositePositionSources_assnlist-class<br>(zzz-DTTCompatibility-class),<br>14                    |
| compositeGroup_assnlist,ArrayDesign_package_methods<br>(zzz-DTTCompatibility-class),<br>14  | CompositeSequence<br>(zzz-DTTCompatibility-class),<br>14  |
| CompositeGroup_assnlist-class   | compositeSequence,CompositeSequence_assnlist-method   |

- (zzz-DTTCompatibility-class),  
14
- compositeSequence, DesignElement\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- CompositeSequence-class  
(zzz-DTTCompatibility-class),  
14
- compositeSequence\_assnlist  
(zzz-DTTCompatibility-class),  
14
- compositeSequence\_assnlist, DesignElement\_package-method  
(zzz-DTTCompatibility-class),  
14
- CompositeSequence\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- compositeSequence\_assnref  
(zzz-DTTCompatibility-class),  
14
- compositeSequence\_assnref, CompositeCompositeMap-method  
(zzz-DTTCompatibility-class),  
14
- compositeSequence\_assnref, ReporterCompositeMap-method  
(zzz-DTTCompatibility-class),  
14
- CompositeSequence\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- compositeSequence\_ref  
(zzz-DTTCompatibility-class),  
14
- compositeSequence\_ref, Composite\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- compositeSequence\_ref, CompositeSequence\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- compositeSequence\_ref, CompositeSequences\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- compositeSequence\_ref, DesignElement\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- CompositeSequence\_ref-class  
(zzz-DTTCompatibility-class),  
14
- compositeSequenceDimension  
(zzz-DTTCompatibility-class),  
14
- compositeSequenceDimension, DesignElementDimension\_assnlist  
(zzz-DTTCompatibility-class),  
14
- CompositeSequenceDimension-class  
(zzz-DTTCompatibility-class),  
14
- compositeSequenceDimension\_ref  
(zzz-DTTCompatibility-class),  
14
- compositeSequenceDimension\_ref, DesignElementDimension\_assnlist  
(zzz-DTTCompatibility-class),  
14
- CompositeSequenceDimension\_ref-class  
(zzz-DTTCompatibility-class),  
14
- compositeSequences\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- compositeSequences\_assnreflist, CompositeGroup-method  
(zzz-DTTCompatibility-class),  
14
- compositeSequences\_assnreflist, CompositeSequenceDimension  
(zzz-DTTCompatibility-class),  
14
- CompositeSequences\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- compound (zzz-DTTCompatibility-class),  
14
- compound, Compound\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Compound-class  
(zzz-DTTCompatibility-class),  
14
- compound\_assnlist  
(zzz-DTTCompatibility-class),  
14
- compound\_assnlist, BioMaterial\_package-method  
(zzz-DTTCompatibility-class),  
14
- Compound\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- compound\_assnref  
(zzz-DTTCompatibility-class),  
14

- 14  
 compound\_assnref,CompoundMeasurement-method  
 (zzz-DTTCompatibility-class),  
 14
- Compound\_assnref-class  
 (zzz-DTTCompatibility-class),  
 14
- compound\_ref  
 (zzz-DTTCompatibility-class),  
 14
- compound\_ref,Compound\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14
- compound\_ref,Labels\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- Compound\_ref-class  
 (zzz-DTTCompatibility-class),  
 14
- compoundIndices\_assnlist  
 (zzz-DTTCompatibility-class),  
 14
- compoundIndices\_assnlist,Compound-method  
 (zzz-DTTCompatibility-class),  
 14
- CompoundIndices\_assnlist-class  
 (zzz-DTTCompatibility-class),  
 14
- compoundMeasurement  
 (zzz-DTTCompatibility-class),  
 14
- compoundMeasurement,ComponentCompounds\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- compoundMeasurement,CompoundMeasurements\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- CompoundMeasurement-class  
 (zzz-DTTCompatibility-class),  
 14
- compoundMeasurements\_assnlist  
 (zzz-DTTCompatibility-class),  
 14
- compoundMeasurements\_assnlist,Treatment-method  
 (zzz-DTTCompatibility-class),  
 14
- CompoundMeasurements\_assnlist-class  
 (zzz-DTTCompatibility-class),
- 14  
 concentrationUnit  
 (zzz-DTTCompatibility-class),  
 14
- concentrationUnit,Unit\_assn-method  
 (zzz-DTTCompatibility-class),  
 14
- ConcentrationUnit-class  
 (zzz-DTTCompatibility-class),  
 14
- ConfidenceIndicator\_assnlist-class  
 (zzz-DTTCompatibility-class),  
 14
- ConfidenceIndicator\_ref-class  
 (zzz-DTTCompatibility-class),  
 14
- confidenceIndicators\_assnreflist  
 (zzz-DTTCompatibility-class),  
 14
- confidenceIndicators\_assnreflist,DerivedSignal-method  
 (zzz-DTTCompatibility-class),  
 14
- confidenceIndicators\_assnreflist,Error-method  
 (zzz-DTTCompatibility-class),  
 14
- confidenceIndicators\_assnreflist,ExpectedValue-method  
 (zzz-DTTCompatibility-class),  
 14
- confidenceIndicators\_assnreflist,Failed-method  
 (zzz-DTTCompatibility-class),  
 14
- confidenceIndicators\_assnreflist,MeasuredSignal-method  
 (zzz-DTTCompatibility-class),  
 14
- confidenceIndicators\_assnreflist,PresentAbsent-method  
 (zzz-DTTCompatibility-class),  
 14
- confidenceIndicators\_assnreflist,PValue-method  
 (zzz-DTTCompatibility-class),  
 14
- confidenceIndicators\_assnreflist,Ratio-method  
 (zzz-DTTCompatibility-class),  
 14
- confidenceIndicators\_assnreflist,SpecializedQuantitationType  
 (zzz-DTTCompatibility-class),  
 14
- ConfidenceIndicators\_assnreflist-class  
 (zzz-DTTCompatibility-class),

- 14
- contact\_assnlist
  - (zzz-DTTCompatibility-class), 14
- contact\_assnlist, AuditAndSecurity\_package-method
  - (zzz-DTTCompatibility-class), 14
- Contact\_assnlist-class
  - (zzz-DTTCompatibility-class), 14
- Contact\_ref-class
  - (zzz-DTTCompatibility-class), 14
- contacts\_assnreflist
  - (zzz-DTTCompatibility-class), 14
- contacts\_assnreflist, Database-method
  - (zzz-DTTCompatibility-class), 14
- Contacts\_assnreflist-class
  - (zzz-DTTCompatibility-class), 14
- containedFeatures\_assnreflist
  - (zzz-DTTCompatibility-class), 14
- containedFeatures\_assnreflist, FeatureDimension-method
  - (zzz-DTTCompatibility-class), 14
- ContainedFeatures\_assnreflist-class
  - (zzz-DTTCompatibility-class), 14
- control (zzz-ARRCompatibility-class), 13
- control, UserAttribute-method
  - (zzz-ARRCompatibility-class), 13
- Control-class
  - (zzz-ARRCompatibility-class), 13
- controlFeatures\_assnreflist
  - (zzz-DTTCompatibility-class), 14
- controlFeatures\_assnreflist, Feature-method
  - (zzz-DTTCompatibility-class), 14
- ControlFeatures\_assnreflist-class
  - (zzz-DTTCompatibility-class), 14
- controlledFeatures\_assnreflist
  - (zzz-DTTCompatibility-class), 14
- controlledFeatures\_assnreflist, Feature-method
  - (zzz-DTTCompatibility-class), 14
- ControlledFeatures\_assnreflist-class
  - (zzz-DTTCompatibility-class), 14
- controlType\_assn
  - (zzz-DTTCompatibility-class), 14
- controlType\_assn, CompositeSequence-method
  - (zzz-DTTCompatibility-class), 14
- controlType\_assn, Feature-method
  - (zzz-DTTCompatibility-class), 14
- controlType\_assn, Reporter-method
  - (zzz-DTTCompatibility-class), 14
- ControlType\_assn-class
  - (zzz-DTTCompatibility-class), 14
- coordinate\_assn
  - (zzz-DTTCompatibility-class), 14
- coordinate\_assn, SeqFeatureLocation-method
  - (zzz-DTTCompatibility-class), 14
- Coordinate\_assn-class
  - (zzz-DTTCompatibility-class), 14
- createdBy (zzz-ARRCompatibility-class), 13
- createdBy, ArraySetFile-method
  - (zzz-ARRCompatibility-class), 13
- createdBy, PhysicalArray-method
  - (zzz-ARRCompatibility-class), 13
- createdBy, TemplateFile-method
  - (zzz-ARRCompatibility-class), 13
- createdDateTime
  - (zzz-ARRCompatibility-class), 13
- createdDateTime, ArraySetFile-method
  - (zzz-ARRCompatibility-class), 13

- 13
- createdDateTime,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- createdDateTime,TemplateFile-method  
(zzz-ARRCompatibility-class),  
13
- createdStep  
(zzz-ARRCompatibility-class),  
13
- createdStep,ArraySetFile-method  
(zzz-ARRCompatibility-class),  
13
- createdStep,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- database (zzz-DTTCompatibility-class),  
14
- database,Database\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Database-class  
(zzz-DTTCompatibility-class),  
14
- database\_assnlist  
(zzz-DTTCompatibility-class),  
14
- database\_assnlist,Description\_package-method  
(zzz-DTTCompatibility-class),  
14
- Database\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- database\_assnref  
(zzz-DTTCompatibility-class),  
14
- database\_assnref,DatabaseEntry-method  
(zzz-DTTCompatibility-class),  
14
- Database\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- database\_ref  
(zzz-DTTCompatibility-class),  
14
- database\_ref,Database\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- Database\_ref-class  
(zzz-DTTCompatibility-class),  
14
- databaseEntry  
(zzz-DTTCompatibility-class),  
14
- databaseEntry,Accessions\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- databaseEntry,DatabaseReferences\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- databaseEntry,ExternalLIMS\_assn-method  
(zzz-DTTCompatibility-class),  
14
- databaseEntry,IdentifierLIMS\_assn-method  
(zzz-DTTCompatibility-class),  
14
- databaseEntry,OntologyReference\_assn-method  
(zzz-DTTCompatibility-class),  
14
- databaseEntry,SequenceDatabases\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- DatabaseEntry-class  
(zzz-DTTCompatibility-class),  
14
- databaseReferences\_assnlist  
(zzz-DTTCompatibility-class),  
14
- databaseReferences\_assnlist,Description-method  
(zzz-DTTCompatibility-class),  
14
- DatabaseReferences\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- dataExternal  
(zzz-DTTCompatibility-class),  
14
- dataExternal,DataExternal\_assn-method  
(zzz-DTTCompatibility-class),  
14
- DataExternal-class  
(zzz-DTTCompatibility-class),  
14
- dataExternal\_assn  
(zzz-DTTCompatibility-class),  
14

- dataExternal\_assn, BioDataCube-method  
(zzz-DTTCompatibility-class),  
[14](#)
- DataExternal\_assn-class  
(zzz-DTTCompatibility-class),  
[14](#)
- dataFormat  
(zzz-DTTCompatibility-class),  
[14](#)
- dataFormat, DataExternal-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataFormatInfoURI  
(zzz-DTTCompatibility-class),  
[14](#)
- dataFormatInfoURI, DataExternal-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataInternal  
(zzz-DTTCompatibility-class),  
[14](#)
- dataInternal, DataInternal\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- DataInternal-class  
(zzz-DTTCompatibility-class),  
[14](#)
- dataInternal\_assn  
(zzz-DTTCompatibility-class),  
[14](#)
- dataInternal\_assn, BioDataCube-method  
(zzz-DTTCompatibility-class),  
[14](#)
- DataInternal\_assn-class  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, DerivedSignal-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, Error-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, ExpectedValue-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, Failed-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, MeasuredSignal-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, NodeValue-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, Parameter-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, PresentAbsent-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, PValue-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, Ratio-method  
(zzz-DTTCompatibility-class),  
[14](#)
- dataType\_assn, SpecializedQuantitationType-method  
(zzz-DTTCompatibility-class),  
[14](#)
- DataType\_assn-class  
(zzz-DTTCompatibility-class),  
[14](#)
- datum (zzz-DTTCompatibility-class), [14](#)
- datum, Datum\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- Datum-class  
(zzz-DTTCompatibility-class),  
[14](#)
- datum\_assn  
(zzz-DTTCompatibility-class),  
[14](#)
- datum\_assn, QuantitationTypeTuple-method  
(zzz-DTTCompatibility-class),  
[14](#)
- Datum\_assn-class  
(zzz-DTTCompatibility-class),  
[14](#)
- defaultValue  
(zzz-ARRCompatibility-class),  
[13](#)
- defaultValue, UserAttribute-method  
(zzz-ARRCompatibility-class),

- 13
- defaultValue\_assn  
(zzz-DTTCompatibility-class),  
14
- defaultValue\_assn,Parameter-method  
(zzz-DTTCompatibility-class),  
14
- DefaultValue\_assn-class  
(zzz-DTTCompatibility-class),  
14
- defectType\_assn  
(zzz-DTTCompatibility-class),  
14
- defectType\_assn,FeatureDefect-method  
(zzz-DTTCompatibility-class),  
14
- defectType\_assn,ZoneDefect-method  
(zzz-DTTCompatibility-class),  
14
- DefectType\_assn-class  
(zzz-DTTCompatibility-class),  
14
- deltaX(zzz-DTTCompatibility-class), 14
- deltaX,PositionDelta-method  
(zzz-DTTCompatibility-class),  
14
- deltaY(zzz-DTTCompatibility-class), 14
- deltaY,PositionDelta-method  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssay  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssay,BioAssay\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssay,DerivedBioAssay\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- DerivedBioAssay-class  
(zzz-DTTCompatibility-class),  
14
- DerivedBioAssay\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssay\_ref  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssay\_ref,BioAssay\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssay\_ref,BioAssayMapTarget\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssay\_ref,BioAssays\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssay\_ref,SourceBioAssays\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssay\_ref,TopLevelBioAssays\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- DerivedBioAssay\_ref-class  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssayData  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssayData,BioAssayData\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssayData,DerivedBioAssayData\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- DerivedBioAssayData-class  
(zzz-DTTCompatibility-class),  
14
- DerivedBioAssayData\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssayData\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssayData\_assnreflist,DerivedBioAssay-method  
(zzz-DTTCompatibility-class),  
14
- DerivedBioAssayData\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssayData\_ref  
(zzz-DTTCompatibility-class),  
14
- derivedBioAssayData\_ref,BioAssayData\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14



|  |  |
|--|--|
| derivedBioAssayData_ref,BioAssayDataSources_assnref-list-method<br>(zzz-DTTCompatibility-class),<br>14       | derivedSignal_ref,TargetQuantitationType_assnref-method<br>(zzz-DTTCompatibility-class),<br>14 |
| derivedBioAssayData_ref,ClusterBioAssayData_assnref-class<br>(zzz-DTTCompatibility-class),<br>14             | derivedSignal_ref-class<br>(zzz-DTTCompatibility-class),<br>14                                 |
| derivedBioAssayData_ref,DerivedBioAssayData_assnref-list-method<br>(zzz-DTTCompatibility-class),<br>14       | description<br>(zzz-DTTCompatibility-class),<br>14   |
| derivedBioAssayData_ref,DerivedBioAssayDataTarget_assnref-list-method<br>(zzz-DTTCompatibility-class),<br>14 | description,DescriptionOptions_assnlist-method<br>(zzz-DTTCompatibility-class),<br>14          |
| DerivedBioAssayData_ref-class<br>(zzz-DTTCompatibility-class),<br>14   | description,NormalizationDescription_assn-method<br>(zzz-DTTCompatibility-class),<br>14        |
| DerivedBioAssayDataTarget_assnref-class<br>(zzz-DTTCompatibility-class),<br>14                               | description,OntologyEntry-method<br>(zzz-DTTCompatibility-class),<br>14                        |
| derivedBioAssayMap_assnreflist<br>(zzz-DTTCompatibility-class),<br>14  | description,QualityControlDescription_assn-method<br>(zzz-DTTCompatibility-class),<br>14       |
| derivedBioAssayMap_assnreflist,DerivedBioAssayMap_assnref-list-method<br>(zzz-DTTCompatibility-class),<br>14 | description,ReplicateDescription_assn-method<br>(zzz-DTTCompatibility-class),<br>14            |
| DerivedBioAssayMap_assnreflist-class<br>(zzz-DTTCompatibility-class),<br>14                                  | Description-class<br>(zzz-DTTCompatibility-class),<br>14                                       |
| derivedSignal<br>(zzz-DTTCompatibility-class),<br>14   | description_package<br>(zzz-DTTCompatibility-class),<br>14                                     |
| derivedSignal,QuantitationType_assnlist-method<br>(zzz-DTTCompatibility-class),<br>14                        | description_package,MAGE-ML-method<br>(zzz-DTTCompatibility-class),<br>14                      |
| DerivedSignal-class<br>(zzz-DTTCompatibility-class),<br>14   | Description_package-class<br>(zzz-DTTCompatibility-class),<br>14                               |
| derivedSignal_ref<br>(zzz-DTTCompatibility-class),<br>14   | descriptions_assnlist<br>(zzz-DTTCompatibility-class),<br>14                                   |
| derivedSignal_ref,QuantitationType_assnref-method<br>(zzz-DTTCompatibility-class),<br>14                     | descriptions_assnlist,Array-method<br>(zzz-DTTCompatibility-class),<br>14                      |
| derivedSignal_ref,QuantitationTypes_assnref-list-method<br>(zzz-DTTCompatibility-class),<br>14               | descriptions_assnlist,ArrayDesign-method<br>(zzz-DTTCompatibility-class),<br>14                |
| derivedSignal_ref,SourcesQuantitationType_assnref-list-method<br>(zzz-DTTCompatibility-class),<br>14         | descriptions_assnlist,ArrayGroup-method<br>(zzz-DTTCompatibility-class),<br>14                 |

descriptions\_assnlist,ArrayManufacture-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,Audit-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,BibliographicReference-description  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,BioAssayCreation-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,BioAssayDataCluster-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,BioAssayDimension-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,BioAssayMap-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,BioAssayTreatment-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,BioSample-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,BioSequence-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,BioSource-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,Channel-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,CompositeCompositeMap-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,CompositeGroup-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,CompositeSequence-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,CompositeSequenceDimension-description  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,Compound-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,Database-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,DerivedBioAssay-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,DerivedBioAssayData-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,DerivedSignal-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,Description-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,Error-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,ExpectedValue-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,Experiment-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,ExperimentalFactor-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,ExperimentDesign-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,FactorValue-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,Failed-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,Feature-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,FeatureDimension-method  
 (zzz-DTTCompatibility-class), 14

descriptions\_assnlist,FeatureExtraction-method  
 (zzz-DTTCompatibility-class), 14

- descriptions\_assnlist,FeatureGroup-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,FeatureReporterMap-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Fiducial-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Hardware-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,HardwareApplication-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Hybridization-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Image-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,ImageAcquisition-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,LabeledExtract-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,MAGE-ML-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,ManufactureLIMS-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,ManufactureLIMSBiomaterial-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,MeasuredBioAssay-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,MeasuredBioAssayData-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,MeasuredSignal-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Node-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,NodeContents-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Organization-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Parameter-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Person-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,PhysicalArrayDesign-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,PhysicalBioAssay-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,PresentAbsent-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Protocol-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,ProtocolApplication-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,PValue-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,QuantitationTypeDimension-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,QuantitationTypeMap-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Ratio-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Reporter-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,ReporterCompositeMap-method  
 (zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,ReporterDimension-method  
 (zzz-DTTCompatibility-class),  
[14](#)

- descriptions\_assnlist,ReporterGroup-method  
(zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Security-method  
(zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,SecurityGroup-method  
(zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,SeqFeature-method  
(zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Software-method  
(zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,SoftwareApplication-method  
(zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,SpecializedQuantitationType-method  
(zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Transformation-method  
(zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Treatment-method  
(zzz-DTTCompatibility-class),  
[14](#)
- descriptions\_assnlist,Zone-method  
(zzz-DTTCompatibility-class),  
[14](#)
- Descriptions\_assnlist-class  
(zzz-DTTCompatibility-class),  
[14](#)
- DesignElement\_assnlist-class  
(zzz-DTTCompatibility-class),  
[14](#)
- designElement\_assnref  
(zzz-DTTCompatibility-class),  
[14](#)
- designElement\_assnref,DesignElementTuple-method  
(zzz-DTTCompatibility-class),  
[14](#)
- DesignElement\_assnref-class  
(zzz-DTTCompatibility-class),  
[14](#)
- designElement\_package  
(zzz-DTTCompatibility-class),  
[14](#)
- designElement\_package,MAGE-ML-method  
(zzz-DTTCompatibility-class),  
[14](#)
- DesignElement\_package-class  
(zzz-DTTCompatibility-class),  
[14](#)
- DesignElement\_ref-class  
(zzz-DTTCompatibility-class),  
[14](#)
- designElementDimension\_assnlist  
(zzz-DTTCompatibility-class),  
[14](#)
- designElementDimension\_assnlist,BioAssayData\_package-method  
(zzz-DTTCompatibility-class),  
[14](#)
- DesignElementDimension\_assnlist-class  
(zzz-DTTCompatibility-class),  
[14](#)
- DesignElementDimension\_assnref  
(zzz-DTTCompatibility-class),  
[14](#)
- designElementDimension\_assnref,DerivedBioAssayData-method  
(zzz-DTTCompatibility-class),  
[14](#)
- designElementDimension\_assnref,MeasuredBioAssayData-method  
(zzz-DTTCompatibility-class),  
[14](#)
- designElementDimension\_assnref,NodeContents-method  
(zzz-DTTCompatibility-class),  
[14](#)
- DesignElementDimension\_assnref-class  
(zzz-DTTCompatibility-class),  
[14](#)
- DesignElementDimension\_ref-class  
(zzz-DTTCompatibility-class),  
[14](#)
- DesignElementMap\_assnlist-class  
(zzz-DTTCompatibility-class),  
[14](#)
- DesignElementMap\_ref-class  
(zzz-DTTCompatibility-class),  
[14](#)
- designElementMapping  
(zzz-DTTCompatibility-class),  
[14](#)
- designElementMapping,DesignElementMapping\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)

- DesignElementMapping-class  
(zzz-DTTCompatibility-class),  
14
- designElementMapping\_assn  
(zzz-DTTCompatibility-class),  
14
- designElementMapping\_assn,Transformation-method  
(zzz-DTTCompatibility-class),  
14
- DesignElementMapping\_assn-class  
(zzz-DTTCompatibility-class),  
14
- designElementMaps\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- designElementMaps\_assnreflist,DesignElementMapping-method  
(zzz-DTTCompatibility-class),  
14
- DesignElementMaps\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- designElementTuple  
(zzz-DTTCompatibility-class),  
14
- designElementTuple,DesignElementTuples\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- DesignElementTuple-class  
(zzz-DTTCompatibility-class),  
14
- designElementTuples\_assnlist  
(zzz-DTTCompatibility-class),  
14
- designElementTuples\_assnlist,BioAssayTuple-method  
(zzz-DTTCompatibility-class),  
14
- DesignElementTuples\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- designProviders\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- designProviders\_assnreflist,ArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- designProviders\_assnreflist,PhysicalArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- DesignProviders\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- directory (accessors), 2
- directory,NetAffxResource-method  
(NetAffxResource-class), 6
- DistanceUnit  
(zzz-DTTCompatibility-class),  
14
- distanceUnit,DistanceUnit\_assn-method  
(zzz-DTTCompatibility-class),  
14
- distanceUnit,Unit\_assn-method  
(zzz-DTTCompatibility-class),  
14
- DistanceUnit-class  
(zzz-DTTCompatibility-class),  
14
- distanceUnit\_assn  
(zzz-DTTCompatibility-class),  
14
- distanceUnit\_assn,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- distanceUnit\_assn,FeatureGroup-method  
(zzz-DTTCompatibility-class),  
14
- distanceUnit\_assn,Fiducial-method  
(zzz-DTTCompatibility-class),  
14
- distanceUnit\_assn,Position-method  
(zzz-DTTCompatibility-class),  
14
- distanceUnit\_assn,PositionDelta-method  
(zzz-DTTCompatibility-class),  
14
- distanceUnit\_assn,Zone-method  
(zzz-DTTCompatibility-class),  
14
- distanceUnit\_assn,ZoneGroup-method  
(zzz-DTTCompatibility-class),  
14
- distanceUnit\_assn,ZoneLayout-method  
(zzz-DTTCompatibility-class),  
14
- DistanceUnit\_assn-class  
(zzz-DTTCompatibility-class),  
14

- DTTCompatibility, 15
- DTTCompatibility-class  
(AffyCompatible-class), 3
- editor (zzz-DTTCompatibility-class), 14
- editor, BibliographicReference-method  
(zzz-DTTCompatibility-class), 14
- email (zzz-DTTCompatibility-class), 14
- email, Organization-method  
(zzz-DTTCompatibility-class), 14
- email, Person-method  
(zzz-DTTCompatibility-class), 14
- error (zzz-DTTCompatibility-class), 14
- error, ConfidenceIndicator\_assnlist-method  
(zzz-DTTCompatibility-class), 14
- error, QuantitationType\_assnlist-method  
(zzz-DTTCompatibility-class), 14
- Error-class  
(zzz-DTTCompatibility-class), 14
- error\_ref (zzz-DTTCompatibility-class), 14
- error\_ref, ConfidenceIndicators\_assnreflist-method  
(zzz-DTTCompatibility-class), 14
- error\_ref, QuantitationType\_assnref-method  
(zzz-DTTCompatibility-class), 14
- error\_ref, QuantitationTypes\_assnreflist-method  
(zzz-DTTCompatibility-class), 14
- error\_ref, SourcesQuantitationType\_assnreflist-method  
(zzz-DTTCompatibility-class), 14
- error\_ref, TargetQuantitationType\_assnref-method  
(zzz-DTTCompatibility-class), 14
- Error\_ref-class  
(zzz-DTTCompatibility-class), 14
- expectedValue  
(zzz-DTTCompatibility-class), 14
- expectedValue, ConfidenceIndicator\_assnlist-method  
(zzz-DTTCompatibility-class), 14
- expectedValue, QuantitationType\_assnlist-method  
(zzz-DTTCompatibility-class), 14
- ExpectedValue-class  
(zzz-DTTCompatibility-class), 14
- expectedValue\_ref  
(zzz-DTTCompatibility-class), 14
- expectedValue\_ref, ConfidenceIndicators\_assnreflist-method  
(zzz-DTTCompatibility-class), 14
- expectedValue\_ref, QuantitationType\_assnref-method  
(zzz-DTTCompatibility-class), 14
- expectedValue\_ref, QuantitationTypes\_assnreflist-method  
(zzz-DTTCompatibility-class), 14
- expectedValue\_ref, SourcesQuantitationType\_assnreflist-method  
(zzz-DTTCompatibility-class), 14
- expectedValue\_ref, TargetQuantitationType\_assnref-method  
(zzz-DTTCompatibility-class), 14
- ExpectedValue\_ref-class  
(zzz-DTTCompatibility-class), 14
- experiment  
(zzz-DTTCompatibility-class), 14
- experiment, Experiment\_assnlist-method  
(zzz-DTTCompatibility-class), 14
- Experiment-class  
(zzz-DTTCompatibility-class), 14
- experiment\_assnlist  
(zzz-DTTCompatibility-class), 14
- experiment\_assnlist, Experiment\_package-method  
(zzz-DTTCompatibility-class), 14
- Experiment\_assnlist-class  
(zzz-DTTCompatibility-class), 14
- experiment\_package

- (zzz-DTTCompatibility-class),  
14
- experiment\_package, MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- Experiment\_package-class  
(zzz-DTTCompatibility-class),  
14
- experimentalFactor  
(zzz-DTTCompatibility-class),  
14
- experimentalFactor, ExperimentalFactors\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ExperimentalFactor-class  
(zzz-DTTCompatibility-class),  
14
- ExperimentalFactor\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- experimentalFactor\_ref  
(zzz-DTTCompatibility-class),  
14
- experimentalFactor\_ref, ExperimentalFactor\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- ExperimentalFactor\_ref-class  
(zzz-DTTCompatibility-class),  
14
- experimentalFactors\_assnlist  
(zzz-DTTCompatibility-class),  
14
- experimentalFactors\_assnlist, ExperimentDesign-method  
(zzz-DTTCompatibility-class),  
14
- ExperimentalFactors\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- experimentDesign  
(zzz-DTTCompatibility-class),  
14
- experimentDesign, ExperimentDesigns\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ExperimentDesign-class  
(zzz-DTTCompatibility-class),  
14
- experimentDesigns\_assnlist  
(zzz-DTTCompatibility-class),  
14
- experimentDesigns\_assnlist, Experiment-method  
(zzz-DTTCompatibility-class),  
14
- ExperimentDesigns\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- exportedFromDB  
(zzz-DTTCompatibility-class),  
14
- exportedFromDB, ExternalReference-method  
(zzz-DTTCompatibility-class),  
14
- exportedFromServer  
(zzz-DTTCompatibility-class),  
14
- exportedFromServer, ExternalReference-method  
(zzz-DTTCompatibility-class),  
14
- exportID (zzz-DTTCompatibility-class),  
14
- exportID, ExternalReference-method  
(zzz-DTTCompatibility-class),  
14
- exportName  
(zzz-DTTCompatibility-class),  
14
- exportName, ExternalReference-method  
(zzz-DTTCompatibility-class),  
14
- externalLIMS\_assn  
(zzz-DTTCompatibility-class),  
14
- externalLIMS\_assn, Compound-method  
(zzz-DTTCompatibility-class),  
14
- ExternalLIMS\_assn-class  
(zzz-DTTCompatibility-class),  
14
- externalReference  
(zzz-DTTCompatibility-class),  
14
- externalReference, ExternalReference\_assn-method  
(zzz-DTTCompatibility-class),  
14
- externalReference\_assn  
(zzz-DTTCompatibility-class),

- 14
- externalReference\_assn,Description-method  
(zzz-DTTCompatibility-class),  
14
- ExternalReference\_assn-class  
(zzz-DTTCompatibility-class),  
14
- factorValue  
(zzz-DTTCompatibility-class),  
14
- factorValue,FactorValues\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- FactorValue-class  
(zzz-DTTCompatibility-class),  
14
- factorValue\_ref  
(zzz-DTTCompatibility-class),  
14
- factorValue\_ref,BioAssayFactorValues\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- FactorValue\_ref-class  
(zzz-DTTCompatibility-class),  
14
- factorValues\_assnlist  
(zzz-DTTCompatibility-class),  
14
- factorValues\_assnlist,ExperimentalFactor-method  
(zzz-DTTCompatibility-class),  
14
- FactorValues\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- failed(zzz-DTTCompatibility-class), 14
- failed,QuantitationType\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Failed-class  
(zzz-DTTCompatibility-class),  
14
- failed\_ref  
(zzz-DTTCompatibility-class),  
14
- failed\_ref,QuantitationType\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- failed\_ref,QuantitationTypes\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- failed\_ref,SourcesQuantitationType\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- failed\_ref,TargetQuantitationType\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- Failed\_ref-class  
(zzz-DTTCompatibility-class),  
14
- failTypes\_assnlist  
(zzz-DTTCompatibility-class),  
14
- failTypes\_assnlist,Reporter-method  
(zzz-DTTCompatibility-class),  
14
- FailTypes\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- fax(zzz-DTTCompatibility-class), 14
- fax,Organization-method  
(zzz-DTTCompatibility-class),  
14
- fax,Person-method  
(zzz-DTTCompatibility-class),  
14
- feature(zzz-DTTCompatibility-class), 14
- feature,DesignElement\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- feature,Features\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Feature-class  
(zzz-DTTCompatibility-class),  
14
- feature\_assnref  
(zzz-DTTCompatibility-class),  
14
- feature\_assnref,FeatureDefect-method  
(zzz-DTTCompatibility-class),  
14
- feature\_assnref,FeatureInformation-method  
(zzz-DTTCompatibility-class),  
14
- feature\_assnref,ManufactureLIMS-method  
(zzz-DTTCompatibility-class),  
14



- 14  
 feature\_assnref,ManufactureLIMSBiomaterial-method  
 (zzz-DTTCompatibility-class),  
 14
- Feature\_assnref-class  
 (zzz-DTTCompatibility-class),  
 14
- feature\_ref  
 (zzz-DTTCompatibility-class),  
 14
- feature\_ref,ContainedFeatures\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- feature\_ref,ControlFeatures\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- feature\_ref,ControlledFeatures\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- feature\_ref,DesignElement\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14
- feature\_ref,Feature\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14
- Feature\_ref-class  
 (zzz-DTTCompatibility-class),  
 14
- featureDefect  
 (zzz-DTTCompatibility-class),  
 14
- featureDefect,FeatureDefects\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- FeatureDefect-class  
 (zzz-DTTCompatibility-class),  
 14
- featureDefects\_assnlist  
 (zzz-DTTCompatibility-class),  
 14
- featureDefects\_assnlist,ArrayManufactureDevialt  
 (zzz-DTTCompatibility-class),  
 14
- FeatureDefects\_assnlist-class  
 (zzz-DTTCompatibility-class),  
 14
- featureDimension  
 (zzz-DTTCompatibility-class),  
 14
- 14  
 featureDimension,DesignElementDimension\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- FeatureDimension-class  
 (zzz-DTTCompatibility-class),  
 14
- featureDimension\_ref  
 (zzz-DTTCompatibility-class),  
 14
- featureDimension\_ref,DesignElementDimension\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14
- FeatureDimension\_ref-class  
 (zzz-DTTCompatibility-class),  
 14
- featureExtraction  
 (zzz-DTTCompatibility-class),  
 14
- featureExtraction,FeatureExtraction\_assn-method  
 (zzz-DTTCompatibility-class),  
 14
- FeatureExtraction-class  
 (zzz-DTTCompatibility-class),  
 14
- featureExtraction\_assn  
 (zzz-DTTCompatibility-class),  
 14
- featureExtraction\_assn,MeasuredBioAssay-method  
 (zzz-DTTCompatibility-class),  
 14
- FeatureExtraction\_assn-class  
 (zzz-DTTCompatibility-class),  
 14
- featureGroup  
 (zzz-DTTCompatibility-class),  
 14
- featureGroup,FeatureGroups\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- featureGroup-class  
 (zzz-DTTCompatibility-class),  
 14
- FeatureGroup\_assnref-class  
 (zzz-DTTCompatibility-class),  
 14
- featureGroup\_ref  
 (zzz-DTTCompatibility-class),  
 14

- 14  
featureGroup\_ref,FeatureGroup\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- FeatureGroup\_ref-class  
(zzz-DTTCompatibility-class),  
14
- featureGroups\_assnlist  
(zzz-DTTCompatibility-class),  
14
- featureGroups\_assnlist,ArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- featureGroups\_assnlist,PhysicalArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- FeatureGroups\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- featureHeight  
(zzz-DTTCompatibility-class),  
14
- featureHeight,FeatureGroup-method  
(zzz-DTTCompatibility-class),  
14
- featureInformation  
(zzz-DTTCompatibility-class),  
14
- featureInformation,FeatureInformationSources\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- FeatureInformation-class  
(zzz-DTTCompatibility-class),  
14
- featureInformationSources\_assnlist  
(zzz-DTTCompatibility-class),  
14
- featureInformationSources\_assnlist,FeatureReporterMap-method  
(zzz-DTTCompatibility-class),  
14
- FeatureInformationSources\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- featureLength  
(zzz-DTTCompatibility-class),  
14
- featureLength,FeatureGroup-method  
(zzz-DTTCompatibility-class),  
14
- 14  
featureIMSSs\_assnlist  
(zzz-DTTCompatibility-class),  
14
- featureIMSSs\_assnlist,ArrayManufacture-method  
(zzz-DTTCompatibility-class),  
14
- FeatureIMSSs\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- featureLocation  
(zzz-DTTCompatibility-class),  
14
- featureLocation,FeatureLocation\_assn-method  
(zzz-DTTCompatibility-class),  
14
- FeatureLocation-class  
(zzz-DTTCompatibility-class),  
14
- featureLocation\_assn  
(zzz-DTTCompatibility-class),  
14
- featureLocation\_assn,Feature-method  
(zzz-DTTCompatibility-class),  
14
- FeatureLocation\_assn-class  
(zzz-DTTCompatibility-class),  
14
- FeatureReporterMap  
(zzz-DTTCompatibility-class),  
14
- featureReporterMap,DesignElementMap\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- featureReporterMap,FeatureReporterMap\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- FeatureReporterMap-class  
(zzz-DTTCompatibility-class),  
14
- featureReporterMap\_assnlist  
(zzz-DTTCompatibility-class),  
14
- featureReporterMap\_assnlist,DesignElement\_package-method  
(zzz-DTTCompatibility-class),  
14
- FeatureReporterMap\_assnlist-class  
(zzz-DTTCompatibility-class),  
14

- 14
- featureReporterMap\_ref  
(zzz-DTTCompatibility-class),  
14
- featureReporterMap\_ref,DesignElementMaps\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- featureReporterMap\_ref,FeatureReporterMaps\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- FeatureReporterMap\_ref-class  
(zzz-DTTCompatibility-class),  
14
- featureReporterMaps\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- featureReporterMaps\_assnreflist,Reporter-method  
(zzz-DTTCompatibility-class),  
14
- FeatureReporterMaps\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- features\_assnlist  
(zzz-DTTCompatibility-class),  
14
- features\_assnlist,FeatureGroup-method  
(zzz-DTTCompatibility-class),  
14
- Features\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- featureShape\_assn  
(zzz-DTTCompatibility-class),  
14
- featureShape\_assn,FeatureGroup-method  
(zzz-DTTCompatibility-class),  
14
- FeatureShape\_assn-class  
(zzz-DTTCompatibility-class),  
14
- featureWidth  
(zzz-DTTCompatibility-class),  
14
- featureWidth,FeatureGroup-method  
(zzz-DTTCompatibility-class),  
14
- fiducial (zzz-DTTCompatibility-class),  
14
- fiducial,Fiducials\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Fiducial-class  
(zzz-DTTCompatibility-class),  
14
- fiducials\_assnlist  
(zzz-DTTCompatibility-class),  
14
- fiducials\_assnlist,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- Fiducials\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- fiducialType\_assn  
(zzz-DTTCompatibility-class),  
14
- fiducialType\_assn,Fiducial-method  
(zzz-DTTCompatibility-class),  
14
- FiducialType\_assn-class  
(zzz-DTTCompatibility-class),  
14
- filenameURI  
(zzz-DTTCompatibility-class),  
14
- filenameURI,DataExternal-method  
(zzz-DTTCompatibility-class),  
14
- firstName (zzz-DTTCompatibility-class),  
14
- firstName,Person-method  
(zzz-DTTCompatibility-class),  
14
- format\_assn  
(zzz-DTTCompatibility-class),  
14
- format\_assn,Image-method  
(zzz-DTTCompatibility-class),  
14
- Format\_assn-class  
(zzz-DTTCompatibility-class),  
14
- guid (zzz-ARRCompatibility-class), 13
- guid,ArraySetFile-method  
(zzz-ARRCompatibility-class),  
13

- guid,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- guid,TemplateFile-method  
(zzz-ARRCompatibility-class),  
13
- hardware (zzz-DTTCompatibility-class),  
14
- hardware,Hardware\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Hardware-class  
(zzz-DTTCompatibility-class),  
14
- hardware\_assnlist  
(zzz-DTTCompatibility-class),  
14
- hardware\_assnlist,Protocol\_package-method  
(zzz-DTTCompatibility-class),  
14
- Hardware\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- hardware\_assnref  
(zzz-DTTCompatibility-class),  
14
- hardware\_assnref,HardwareApplication-method  
(zzz-DTTCompatibility-class),  
14
- hardware\_assnref,Software-method  
(zzz-DTTCompatibility-class),  
14
- Hardware\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- hardware\_ref  
(zzz-DTTCompatibility-class),  
14
- hardware\_ref,Hardware\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- hardware\_ref,Hardwares\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- Hardware\_ref-class  
(zzz-DTTCompatibility-class),  
14
- hardwareApplication  
(zzz-DTTCompatibility-class),  
14
- hardwareApplication,HardwareApplications\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- HardwareApplication-class  
(zzz-DTTCompatibility-class),  
14
- hardwareApplications\_assnlist  
(zzz-DTTCompatibility-class),  
14
- hardwareApplications\_assnlist,ProtocolApplication-method  
(zzz-DTTCompatibility-class),  
14
- HardwareApplications\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- hardwareManufacturers\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- hardwareManufacturers\_assnreflist,Hardware-method  
(zzz-DTTCompatibility-class),  
14
- HardwareManufacturers\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- hardwares\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- hardwares\_assnreflist,Protocol-method  
(zzz-DTTCompatibility-class),  
14
- Hardwares\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- higherLevelAnalysis\_package  
(zzz-DTTCompatibility-class),  
14
- higherLevelAnalysis\_package,MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- HigherLevelAnalysis\_package-class  
(zzz-DTTCompatibility-class),  
14
- hybridization  
(zzz-DTTCompatibility-class),  
14
- hybridization,BioAssayCreation\_assn-method

- (zzz-DTTCompatibility-class),  
14
- Hybridization-class  
(zzz-DTTCompatibility-class),  
14
- identifier  
(zzz-DTTCompatibility-class),  
14
- identifier,Array-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Array\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ArrayDesign\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ArrayGroup\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ArrayManufacture-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ArrayManufacture\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioAssay\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioAssayCreation-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioAssayData\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioAssayDataCluster-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioAssayDataCluster\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioAssayDimension-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioAssayDimension\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioAssayMap-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioAssayMap\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioAssayTreatment-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioMaterial\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioSample-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioSample\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioSequence-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioSequence\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioSource-method  
(zzz-DTTCompatibility-class),  
14
- identifier,BioSource\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Channel-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Channel\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,CompositeCompositeMap-method  
(zzz-DTTCompatibility-class),  
14
- identifier,CompositeCompositeMap\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,CompositeGroup-method

- (zzz-DTTCompatibility-class),  
14
- identifier, CompositeGroup\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, CompositeSequence-method  
(zzz-DTTCompatibility-class),  
14
- identifier, CompositeSequence\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, CompositeSequenceDimension-method  
(zzz-DTTCompatibility-class),  
14
- identifier, CompositeSequenceDimension\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Compound-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Compound\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, ConfidenceIndicator\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Contact\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Database-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Database\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, DerivedBioAssay-method  
(zzz-DTTCompatibility-class),  
14
- identifier, DerivedBioAssay\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, DerivedBioAssayData-method  
(zzz-DTTCompatibility-class),  
14
- identifier, DerivedBioAssayData\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, DerivedSignal-method  
(zzz-DTTCompatibility-class),  
14
- identifier, DerivedSignal\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, DesignElement\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, DesignElementDimension\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, DesignElementMap\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Error-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Error\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, ExpectedValue-method  
(zzz-DTTCompatibility-class),  
14
- identifier, ExpectedValue\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Experiment-method  
(zzz-DTTCompatibility-class),  
14
- identifier, ExperimentalFactor-method  
(zzz-DTTCompatibility-class),  
14
- identifier, ExperimentalFactor\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, FactorValue-method  
(zzz-DTTCompatibility-class),  
14
- identifier, FactorValue\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Failed-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Failed\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier, Feature-method

- (zzz-DTTCompatibility-class),  
[14](#)
- identifier, Feature\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, FeatureDimension-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, FeatureDimension\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, FeatureExtraction-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, FeatureGroup-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, FeatureGroup\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, FeatureReporterMap-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, FeatureReporterMap\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Hardware-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Hardware\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Hybridization-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Image-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Image\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, ImageAcquisition-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, LabeledExtract-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, LabeledExtract\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, MAGE-ML-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, MeasuredBioAssay-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, MeasuredBioAssay\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, MeasuredBioAssayData-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, MeasuredBioAssayData\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, MeasuredSignal-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, MeasuredSignal\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Organization-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Organization\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Parameter-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Parameter\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Person-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, Person\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, PhysicalArrayDesign-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, PhysicalArrayDesign\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- identifier, PhysicalBioAssay-method

- (zzz-DTTCompatibility-class),  
14
- identifier,PhysicalBioAssay\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,PresentAbsent-method  
(zzz-DTTCompatibility-class),  
14
- identifier,PresentAbsent\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Protocol-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Protocol\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,PValue-method  
(zzz-DTTCompatibility-class),  
14
- identifier,PValue\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,QuantitationType\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,QuantitationTypeDimension-method  
(zzz-DTTCompatibility-class),  
14
- identifier,QuantitationTypeDimension\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,QuantitationTypeMap-method  
(zzz-DTTCompatibility-class),  
14
- identifier,QuantitationTypeMap\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Ratio-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Ratio\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Reporter-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Reporter\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ReporterCompositeMap-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ReporterCompositeMap\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ReporterDimension-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ReporterDimension\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ReporterGroup-method  
(zzz-DTTCompatibility-class),  
14
- identifier,ReporterGroup\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Security-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Security\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,SecurityGroup-method  
(zzz-DTTCompatibility-class),  
14
- identifier,SecurityGroup\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Software-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Software\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,SpecializedQuantitationType-method  
(zzz-DTTCompatibility-class),  
14
- identifier,SpecializedQuantitationType\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,StandardQuantitationType\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Transformation-method  
(zzz-DTTCompatibility-class),  
14



- (zzz-DTTCompatibility-class),  
14
- identifier,Treatment-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Zone-method  
(zzz-DTTCompatibility-class),  
14
- identifier,Zone\_ref-method  
(zzz-DTTCompatibility-class),  
14
- identifierLIMS\_assn  
(zzz-DTTCompatibility-class),  
14
- identifierLIMS\_assn,ManufactureLIMS-method  
(zzz-DTTCompatibility-class),  
14
- identifierLIMS\_assn,ManufactureLIMSBiomaterial-method  
(zzz-DTTCompatibility-class),  
14
- IdentifierLIMS\_assn-class  
(zzz-DTTCompatibility-class),  
14
- Image-class  
(zzz-DTTCompatibility-class),  
14
- image\_ref (zzz-DTTCompatibility-class),  
14
- image\_ref,Images\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- Image\_ref-class  
(zzz-DTTCompatibility-class),  
14
- imageAcquisition  
(zzz-DTTCompatibility-class),  
14
- imageAcquisition,BioAssayTreatments\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ImageAcquisition-class  
(zzz-DTTCompatibility-class),  
14
- images\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- images\_assnreflist,ImageAcquisition-method  
(zzz-DTTCompatibility-class),  
14
- 14
- Images\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- immobilizedCharacteristics\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- immobilizedCharacteristics\_assnreflist,Reporter-method  
(zzz-DTTCompatibility-class),  
14
- ImmobilizedCharacteristics\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- information\_assnref  
(zzz-DTTCompatibility-class),  
14
- information\_assnref,Array-method  
(zzz-DTTCompatibility-class),  
14
- Information\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- Internal-class, 4
- isApproximateLength  
(zzz-DTTCompatibility-class),  
14
- isApproximateLength,BioSequence-method  
(zzz-DTTCompatibility-class),  
14
- isBackground  
(zzz-DTTCompatibility-class),  
14
- isBackground,DerivedSignal-method  
(zzz-DTTCompatibility-class),  
14
- isBackground,Error-method  
(zzz-DTTCompatibility-class),  
14
- isBackground,ExpectedValue-method  
(zzz-DTTCompatibility-class),  
14
- isBackground,Failed-method  
(zzz-DTTCompatibility-class),  
14
- isBackground,MeasuredSignal-method  
(zzz-DTTCompatibility-class),  
14
- isBackground,PresentAbsent-method

- (zzz-DTTCompatibility-class),  
14
- isBackground,PValue-method  
(zzz-DTTCompatibility-class),  
14
- isBackground,Ratio-method  
(zzz-DTTCompatibility-class),  
14
- isBackground,SpecializedQuantitationType-method  
(zzz-DTTCompatibility-class),  
14
- isCircular  
(zzz-DTTCompatibility-class),  
14
- isCircular,BioSequence-method  
(zzz-DTTCompatibility-class),  
14
- isSolvent(zzz-DTTCompatibility-class),  
14
- isSolvent,Compound-method  
(zzz-DTTCompatibility-class),  
14
- issue(zzz-DTTCompatibility-class), 14
- issue,BibliographicReference-method  
(zzz-DTTCompatibility-class),  
14
- kindCV(zzz-DTTCompatibility-class), 14
- kindCV,Measurement-method  
(zzz-DTTCompatibility-class),  
14
- labeledExtract  
(zzz-DTTCompatibility-class),  
14
- labeledExtract,BioMaterial\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- LabeledExtract-class  
(zzz-DTTCompatibility-class),  
14
- labeledExtract\_ref  
(zzz-DTTCompatibility-class),  
14
- labeledExtract\_ref,BioMaterial\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- LabeledExtract\_ref-class  
(zzz-DTTCompatibility-class),
- 14
- labels\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- labels\_assnreflist,Channel-method  
(zzz-DTTCompatibility-class),  
14
- labels\_assnreflist,LabeledExtract-method  
(zzz-DTTCompatibility-class),  
14
- Labels\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- lastName(zzz-DTTCompatibility-class),  
14
- lastName,Person-method  
(zzz-DTTCompatibility-class),  
14
- length,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- length,BioSequence-method  
(zzz-DTTCompatibility-class),  
14
- libraryPackageName  
(zzz-ARRCompatibility-class),  
13
- libraryPackageName,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- list, 4
- lowerRightX  
(zzz-DTTCompatibility-class),  
14
- lowerRightX,Zone-method  
(zzz-DTTCompatibility-class),  
14
- lowerRightY  
(zzz-DTTCompatibility-class),  
14
- lowerRightY,Zone-method  
(zzz-DTTCompatibility-class),  
14
- MAGE-ML-class  
(zzz-DTTCompatibility-class),  
14
- make(zzz-DTTCompatibility-class), 14
- make,Hardware-method

- (zzz-DTTCompatibility-class),  
14
- manufactureLIMS  
(zzz-DTTCompatibility-class),  
14
- manufactureLIMS, FeatureLIMSS\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ManufactureLIMS-class  
(zzz-DTTCompatibility-class),  
14
- manufactureLIMSBiomaterial  
(zzz-DTTCompatibility-class),  
14
- manufactureLIMSBiomaterial, FeatureLIMSS\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ManufactureLIMSBiomaterial-class  
(zzz-DTTCompatibility-class),  
14
- manufacturingDate  
(zzz-DTTCompatibility-class),  
14
- manufacturingDate, ArrayManufacture-method  
(zzz-DTTCompatibility-class),  
14
- massUnit (zzz-DTTCompatibility-class),  
14
- massUnit, Unit\_assn-method  
(zzz-DTTCompatibility-class),  
14
- MassUnit-class  
(zzz-DTTCompatibility-class),  
14
- masterFileGUID  
(zzz-ARRCompatibility-class),  
13
- masterFileGUID, PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- masterFileName  
(zzz-ARRCompatibility-class),  
13
- masterFileName, PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- materialType\_assn  
(zzz-DTTCompatibility-class),  
14
- materialType\_assn, BioSample-method  
(zzz-DTTCompatibility-class),  
14
- materialType\_assn, BioSource-method  
(zzz-DTTCompatibility-class),  
14
- materialType\_assn, LabeledExtract-method  
(zzz-DTTCompatibility-class),  
14
- MaterialType\_assn-class  
(zzz-DTTCompatibility-class),  
14
- measuredBioAssay  
(zzz-DTTCompatibility-class),  
14
- measuredBioAssay, BioAssay\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- measuredBioAssay, MeasuredBioAssay\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- MeasuredBioAssay-class  
(zzz-DTTCompatibility-class),  
14
- MeasuredBioAssay\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- measuredBioAssay\_ref  
(zzz-DTTCompatibility-class),  
14
- measuredBioAssay\_ref, BioAssay\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- measuredBioAssay\_ref, BioAssays\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- measuredBioAssay\_ref, MeasuredBioAssayTarget\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- measuredBioAssay\_ref, SourceBioAssays\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- measuredBioAssay\_ref, TopLevelBioAssays\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- MeasuredBioAssay\_ref-class  
(zzz-DTTCompatibility-class),

- 14  
 measuredBioAssayData  
 (zzz-DTTCompatibility-class),  
 14
- measuredBioAssayData,BioAssayData\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- measuredBioAssayData,MeasuredBioAssayData\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- MeasuredBioAssayData-class  
 (zzz-DTTCompatibility-class),  
 14
- MeasuredBioAssayData\_assnlist-class  
 (zzz-DTTCompatibility-class),  
 14
- measuredBioAssayData\_assnreflist  
 (zzz-DTTCompatibility-class),  
 14
- measuredBioAssayData\_assnreflist,MeasuredBioAssayData\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- MeasuredBioAssayData\_assnreflist-class  
 (zzz-DTTCompatibility-class),  
 14
- measuredBioAssayData\_ref  
 (zzz-DTTCompatibility-class),  
 14
- measuredBioAssayData\_ref,BioAssayData\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- measuredBioAssayData\_ref,BioAssayDataSources\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- measuredBioAssayData\_ref,ClusterBioAssayData\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- measuredBioAssayData\_ref,MeasuredBioAssayData\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- MeasuredBioAssayData\_ref-class  
 (zzz-DTTCompatibility-class),  
 14
- MeasuredBioAssayTarget\_assnref-class  
 (zzz-DTTCompatibility-class),  
 14
- measuredSignal  
 (zzz-DTTCompatibility-class),  
 14
- measuredSignal,QuantitationType\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- measuredSignal-class  
 (zzz-DTTCompatibility-class),  
 14
- measuredSignal\_ref  
 (zzz-DTTCompatibility-class),  
 14
- measuredSignal\_ref,QuantitationType\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14
- measuredSignal\_ref,QuantitationTypes\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- measuredSignal\_ref,SourcesQuantitationType\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- measuredSignal\_ref,TargetQuantitationType\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14
- MeasuredSignal\_ref-class  
 (zzz-DTTCompatibility-class),  
 14
- measurement  
 (zzz-DTTCompatibility-class),  
 14
- measurement,ActionMeasurement\_assn-method  
 (zzz-DTTCompatibility-class),  
 14
- measurement,DefaultValue\_assn-method  
 (zzz-DTTCompatibility-class),  
 14
- measurement,Measurement\_assn-method  
 (zzz-DTTCompatibility-class),  
 14
- Measurement-class  
 (zzz-DTTCompatibility-class),  
 14
- measurement\_assn  
 (zzz-DTTCompatibility-class),  
 14
- measurement\_assn,BioMaterialMeasurement-method  
 (zzz-DTTCompatibility-class),  
 14
- measurement\_assn,CompoundMeasurement-method  
 (zzz-DTTCompatibility-class),  
 14

- 14
- measurement\_assn,FactorValue-method  
(zzz-DTTCompatibility-class),  
14
- Measurement\_assn-class  
(zzz-DTTCompatibility-class),  
14
- measurement\_package  
(zzz-DTTCompatibility-class),  
14
- measurement\_package,MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- Measurement\_package-class  
(zzz-DTTCompatibility-class),  
14
- mediaCol (zzz-ARRCompatibility-class),  
13
- mediaCol,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- mediaFileGUID  
(zzz-ARRCompatibility-class),  
13
- mediaFileGUID,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- mediaFileName  
(zzz-ARRCompatibility-class),  
13
- mediaFileName,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- mediaRow (zzz-ARRCompatibility-class),  
13
- mediaRow,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- mediaType (zzz-ARRCompatibility-class),  
13
- mediaType,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- members\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- members\_assnreflist,SecurityGroup-method  
(zzz-DTTCompatibility-class),  
14
- Members\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- midInitials  
(zzz-DTTCompatibility-class),  
14
- midInitials,Person-method  
(zzz-DTTCompatibility-class),  
14
- mismatchInformation  
(zzz-DTTCompatibility-class),  
14
- mismatchInformation,MismatchInformation\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- MismatchInformation-class  
(zzz-DTTCompatibility-class),  
14
- mismatchInformation\_assnlist  
(zzz-DTTCompatibility-class),  
14
- mismatchInformation\_assnlist,CompositePosition-method  
(zzz-DTTCompatibility-class),  
14
- mismatchInformation\_assnlist,FeatureInformation-method  
(zzz-DTTCompatibility-class),  
14
- mismatchInformation\_assnlist,ReporterPosition-method  
(zzz-DTTCompatibility-class),  
14
- MismatchInformation\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- model (zzz-DTTCompatibility-class), 14
- model,Hardware-method  
(zzz-DTTCompatibility-class),  
14
- name (zzz-DTTCompatibility-class), 14
- name,Array-method  
(zzz-DTTCompatibility-class),  
14
- name,Array\_ref-method  
(zzz-DTTCompatibility-class),  
14
- name,ArrayAttribute-method  
(zzz-ARRCompatibility-class),  
13

name,ArrayDesign-method  
     (zzz-DTTCompatibility-class),  
     14

name,ArrayDesign\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,ArrayGroup-method  
     (zzz-DTTCompatibility-class),  
     14

name,ArrayGroup\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,ArrayManufacture-method  
     (zzz-DTTCompatibility-class),  
     14

name,ArrayManufacture\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioAssay\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioAssayCreation-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioAssayData\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioAssayDataCluster-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioAssayDataCluster\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioAssayDimension-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioAssayDimension\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioAssayMap-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioAssayMap\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioAssayTreatment-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioMaterial\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioSample-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioSample\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioSequence-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioSequence\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioSource-method  
     (zzz-DTTCompatibility-class),  
     14

name,BioSource\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,Channel-method  
     (zzz-DTTCompatibility-class),  
     14

name,Channel\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,CompositeCompositeMap-method  
     (zzz-DTTCompatibility-class),  
     14

name,CompositeCompositeMap\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,CompositeGroup-method  
     (zzz-DTTCompatibility-class),  
     14

name,CompositeGroup\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,CompositeSequence-method  
     (zzz-DTTCompatibility-class),  
     14

name,CompositeSequence\_ref-method  
     (zzz-DTTCompatibility-class),  
     14

name,CompositeSequenceDimension-method  
     (zzz-DTTCompatibility-class),  
     14

name, CompositeSequenceDimension\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Compound-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Compound\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, ConfidenceIndicator\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Contact\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Database-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Database\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, DerivedBioAssay-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, DerivedBioAssay\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, DerivedBioAssayData-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, DerivedBioAssayData\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, DerivedSignal-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, DerivedSignal\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, DesignElement\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, DesignElementDimension\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, DesignElementMap\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Error-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Error\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, ExpectedValue-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, ExpectedValue\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Experiment-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, ExperimentalFactor-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, ExperimentalFactor\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, FactorValue-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, FactorValue\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Failed-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Failed\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Feature-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, Feature\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, FeatureDimension-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, FeatureDimension\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)

name, FeatureExtraction-method  
(zzz-DTTCompatibility-class),  
[14](#)

- name, FeatureGroup-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, FeatureGroup\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, FeatureReporterMap-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, FeatureReporterMap\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Hardware-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Hardware\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Hybridization-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Image-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Image\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, ImageAcquisition-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, LabeledExtract-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, LabeledExtract\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, MAGE-ML-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, MeasuredBioAssay-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, MeasuredBioAssay\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, MeasuredBioAssayData-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, MeasuredBioAssayData\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, MeasuredSignal-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, MeasuredSignal\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, NameValueType-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, NodeValue-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Organization-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Organization\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Parameter-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Parameter\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Person-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Person\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, PhysicalArrayDesign-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, PhysicalArrayDesign\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, PhysicalBioAssay-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, PhysicalBioAssay\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, PresentAbsent-method  
(zzz-DTTCompatibility-class),  
[14](#)



- name, PresentAbsent\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Protocol-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Protocol\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, PValue-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, PValue\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, QuantitationType\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, QuantitationTypeDimension-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, QuantitationTypeDimension\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, QuantitationTypeMap-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, QuantitationTypeMap\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Ratio-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Ratio\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Reporter-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Reporter\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, ReporterCompositeMap-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, ReporterCompositeMap\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, ReporterDimension-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, ReporterDimension\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, ReporterGroup-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, ReporterGroup\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Security-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Security\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, SecurityGroup-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, SecurityGroup\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Software-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Software\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, SpecializedQuantitationType-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, SpecializedQuantitationType\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, StandardQuantitationType\_ref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Transformation-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, Treatment-method  
(zzz-DTTCompatibility-class),  
[14](#)
- name, UserAttribute-method  
(zzz-ARRCompatibility-class),  
[13](#)

name, Zone-method  
     (zzz-DTTCompatibility-class),  
     14  
 name, Zone\_ref-method  
     (zzz-DTTCompatibility-class),  
     14  
 names, AffxArray-method  
     (zzz-NetAffxCompatibility-class),  
     16  
 names, NetAffxResource-method  
     (NetAffxResource-class), 6  
 namespace (zzz-ARRCompatibility-class),  
     13  
 namespace, UserAttribute-method  
     (zzz-ARRCompatibility-class),  
     13  
 nameValueType  
     (zzz-DTTCompatibility-class),  
     14  
 nameValueType, PropertySets\_assnlist-method  
     (zzz-DTTCompatibility-class),  
     14  
 nameValueType, QualityControlStatistics\_assnlist-method  
     (zzz-DTTCompatibility-class),  
     14  
 nameValueType, SummaryStatistics\_assnlist-method  
     (zzz-DTTCompatibility-class),  
     14  
 NameValueType-class  
     (zzz-DTTCompatibility-class),  
     14  
 NetAffxCompatibility, 6, 16  
 NetAffxCompatibility-class  
     (AffyCompatible-class), 3  
 NetAffxResource, 3, 4, 5, 10, 16, 17  
 NetAffxResource-accessors (accessors), 2  
 NetAffxResource-class, 6  
 newSequence  
     (zzz-DTTCompatibility-class),  
     14  
 newSequence, MismatchInformation-method  
     (zzz-DTTCompatibility-class),  
     14  
 node (zzz-DTTCompatibility-class), 14  
 node, Nodes\_assnlist-method  
     (zzz-DTTCompatibility-class),  
     14  
 Node-class  
     (zzz-DTTCompatibility-class),  
     14  
 nodeContents  
     (zzz-DTTCompatibility-class),  
     14  
 nodeContents, NodeContents\_assnlist-method  
     (zzz-DTTCompatibility-class),  
     14  
 NodeContents-class  
     (zzz-DTTCompatibility-class),  
     14  
 nodeContents\_assnlist  
     (zzz-DTTCompatibility-class),  
     14  
 nodeContents\_assnlist, Node-method  
     (zzz-DTTCompatibility-class),  
     14  
 NodeContents\_assnlist-class  
     (zzz-DTTCompatibility-class),  
     14  
 nodes\_assnlist  
     (zzz-DTTCompatibility-class),  
     14  
 nodes\_assnlist, BioAssayDataCluster-method  
     (zzz-DTTCompatibility-class),  
     14  
 nodes\_assnlist, Node-method  
     (zzz-DTTCompatibility-class),  
     14  
 Nodes\_assnlist-class  
     (zzz-DTTCompatibility-class),  
     14  
 nodeValue (zzz-DTTCompatibility-class),  
     14  
 nodeValue, NodeValue\_assnlist-method  
     (zzz-DTTCompatibility-class),  
     14  
 NodeValue-class  
     (zzz-DTTCompatibility-class),  
     14  
 nodeValue\_assnlist  
     (zzz-DTTCompatibility-class),  
     14  
 nodeValue\_assnlist, Node-method  
     (zzz-DTTCompatibility-class),  
     14  
 NodeValue\_assnlist-class  
     (zzz-DTTCompatibility-class),  
     14

- 14
- normalizationDescription\_assn  
(zzz-DTTCompatibility-class),  
14
- normalizationDescription\_assn,ExperimentDesign-method(zzz-DTTCompatibility-class),  
(zzz-DTTCompatibility-class),  
14
- NormalizationDescription\_assn-class  
(zzz-DTTCompatibility-class),  
14
- numArrays (zzz-DTTCompatibility-class),  
14
- numArrays,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- numberOfFeatures  
(zzz-DTTCompatibility-class),  
14
- numberOfFeatures,ArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- numberOfFeatures,PhysicalArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- numFeaturesPerCol  
(zzz-DTTCompatibility-class),  
14
- numFeaturesPerCol,ZoneLayout-method  
(zzz-DTTCompatibility-class),  
14
- numFeaturesPerRow  
(zzz-DTTCompatibility-class),  
14
- numFeaturesPerRow,ZoneLayout-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntries\_assnlist  
(zzz-DTTCompatibility-class),  
14
- ontologyEntries\_assnlist,BioSequence-method  
(zzz-DTTCompatibility-class),  
14
- OntologyEntries\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,Action\_assn-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,Annotations\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,Associations\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,Category\_assn-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,Characteristics\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,CompoundIndices\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,ControlType\_assn-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,DataType\_assn-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,DefectType\_assn-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,FailTypes\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,FeatureShape\_assn-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,FiducialType\_assn-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,Format\_assn-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,MaterialType\_assn-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,OntologyEntries\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ontologyEntry,Parameters\_assnlist-method  
(zzz-DTTCompatibility-class),  
14

- ontologyEntry, PolymerType\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyEntry, Roles\_assnlist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyEntry, Scale\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyEntry, Species\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyEntry, SubstrateType\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyEntry, SurfaceType\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyEntry, TechnologyType\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyEntry, Type\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyEntry, Types\_assnlist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyEntry, Value\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyEntry, WarningType\_assn-method  
(zzz-DTTCompatibility-class),  
[14](#)
- OntologyEntry-class  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyReference\_assn  
(zzz-DTTCompatibility-class),  
[14](#)
- ontologyReference\_assn, OntologyEntry-method  
(zzz-DTTCompatibility-class),  
[14](#)
- OntologyReference\_assn-class  
(zzz-DTTCompatibility-class),  
[14](#)
- organization  
(zzz-DTTCompatibility-class),  
[14](#)
- organization, Contact\_assnlist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization, Organization\_assnlist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- Organization-class  
(zzz-DTTCompatibility-class),  
[14](#)
- Organization\_assnlist-class  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, Affiliation\_assnref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, ArrayManufacturers\_assnreflist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, Contacts\_assnreflist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, DesignProviders\_assnreflist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, HardwareManufacturers\_assnreflist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, Members\_assnreflist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, Owner\_assnreflist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, Parent\_assnref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, Performer\_assnref-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, Providers\_assnreflist-method  
(zzz-DTTCompatibility-class),  
[14](#)
- organization\_ref, SoftwareManufacturers\_assnreflist-method  
(zzz-DTTCompatibility-class),  
[14](#)

- organization\_ref,SourceContact\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- Organization\_ref-class  
(zzz-DTTCompatibility-class),  
14
- orientationMark  
(zzz-DTTCompatibility-class),  
14
- orientationMark,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- orientationMarkPosition  
(zzz-DTTCompatibility-class),  
14
- orientationMarkPosition,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- originalProjectName  
(zzz-ARRCompatibility-class),  
13
- originalProjectName,ArraySetFile-method  
(zzz-ARRCompatibility-class),  
13
- originRelativeTo  
(zzz-DTTCompatibility-class),  
14
- originRelativeTo,Array-method  
(zzz-DTTCompatibility-class),  
14
- otherKind (zzz-DTTCompatibility-class),  
14
- otherKind,Measurement-method  
(zzz-DTTCompatibility-class),  
14
- owner\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- owner\_assnreflist,Security-method  
(zzz-DTTCompatibility-class),  
14
- Owner\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- pages (zzz-DTTCompatibility-class), 14
- pages,BibliographicReference-method  
(zzz-DTTCompatibility-class),  
14
- parameter (zzz-DTTCompatibility-class),  
14
- parameter,ParameterTypes\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Parameter-class  
(zzz-DTTCompatibility-class),  
14
- parameter\_ref  
(zzz-DTTCompatibility-class),  
14
- parameter\_ref,ParameterType\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- Parameter\_ref-class  
(zzz-DTTCompatibility-class),  
14
- parameters\_assnlist  
(zzz-DTTCompatibility-class),  
14
- parameters\_assnlist,BibliographicReference-method  
(zzz-DTTCompatibility-class),  
14
- Parameters\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- parameterType\_assnref  
(zzz-DTTCompatibility-class),  
14
- parameterType\_assnref,ParameterValue-method  
(zzz-DTTCompatibility-class),  
14
- ParameterType\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- parameterTypes\_assnlist  
(zzz-DTTCompatibility-class),  
14
- parameterTypes\_assnlist,Hardware-method  
(zzz-DTTCompatibility-class),  
14
- parameterTypes\_assnlist,Protocol-method  
(zzz-DTTCompatibility-class),  
14
- parameterTypes\_assnlist,Software-method  
(zzz-DTTCompatibility-class),  
14
- ParameterTypes\_assnlist-class

- (zzz-DTTCompatibility-class),  
14
- parameterValue  
(zzz-DTTCompatibility-class),  
14
- parameterValue,ParameterValues\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ParameterValue-class  
(zzz-DTTCompatibility-class),  
14
- parameterValues\_assnlist  
(zzz-DTTCompatibility-class),  
14
- parameterValues\_assnlist,HardwareApplication-Method  
(zzz-DTTCompatibility-class),  
14
- parameterValues\_assnlist,ProtocolApplication-Method  
(zzz-DTTCompatibility-class),  
14
- parameterValues\_assnlist,SoftwareApplication-Method  
(zzz-DTTCompatibility-class),  
14
- ParameterValues\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- parent\_assnref  
(zzz-DTTCompatibility-class),  
14
- parent\_assnref,Organization-method  
(zzz-DTTCompatibility-class),  
14
- Parent\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- password (accessors), 2
- password,NetAffxResource-method  
(NetAffxResource-class), 6
- patassignmentMethod  
(zzz-ARRCompatibility-class),  
13
- patassignmentMethod,PhysicalArray-method  
(zzz-ARRCompatibility-class),  
13
- performer\_assnref  
(zzz-DTTCompatibility-class),  
14
- performer\_assnref,Audit-method  
(zzz-DTTCompatibility-class),  
14
- (zzz-DTTCompatibility-class),  
14
- Performer\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- performers\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- performers\_assnreflist,ProtocolApplication-method  
(zzz-DTTCompatibility-class),  
14
- Performers\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- Performers\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- person\_ref,ArrayManufacturers\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- person\_ref,Contacts\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- person\_ref,DesignProviders\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- person\_ref,HardwareManufacturers\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- person\_ref,Members\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- person\_ref,Owner\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- person\_ref,Performer\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- person\_ref,Performers\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- person\_ref,Providers\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14

- (zzz-DTTCompatibility-class),  
14
- person\_ref, SoftwareManufacturers\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- person\_ref, SourceContact\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- Person\_ref-class  
(zzz-DTTCompatibility-class),  
14
- phone (zzz-DTTCompatibility-class), 14
- phone, Organization-method  
(zzz-DTTCompatibility-class),  
14
- phone, Person-method  
(zzz-DTTCompatibility-class),  
14
- physicalArray  
(zzz-ARRCompatibility-class),  
13
- physicalArray, PhysicalArrays-method  
(zzz-ARRCompatibility-class),  
13
- PhysicalArray-class  
(zzz-ARRCompatibility-class),  
13
- physicalArrayDesign  
(zzz-DTTCompatibility-class),  
14
- physicalArrayDesign, ArrayDesign\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- PhysicalArrayDesign-class  
(zzz-DTTCompatibility-class),  
14
- physicalArrayDesign\_ref  
(zzz-DTTCompatibility-class),  
14
- physicalArrayDesign\_ref, ArrayDesign\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- PhysicalArrayDesign\_ref-class  
(zzz-DTTCompatibility-class),  
14
- physicalArrays  
(zzz-ARRCompatibility-class),  
13
- physicalArrays, ArraySetFile-method  
(zzz-ARRCompatibility-class),  
13
- PhysicalArrays-class  
(zzz-ARRCompatibility-class),  
13
- physicalBioAssay  
(zzz-DTTCompatibility-class),  
14
- physicalBioAssay, BioAssay\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- physicalBioAssay, PhysicalBioAssay\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- PhysicalBioAssay-class  
(zzz-DTTCompatibility-class),  
14
- PhysicalBioAssay\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- PhysicalBioAssay\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- physicalBioAssay\_ref  
(zzz-DTTCompatibility-class),  
14
- physicalBioAssay\_ref, BioAssay\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- physicalBioAssay\_ref, BioAssays\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- physicalBioAssay\_ref, PhysicalBioAssay\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- physicalBioAssay\_ref, PhysicalBioAssaySource\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- physicalBioAssay\_ref, PhysicalBioAssayTarget\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- physicalBioAssay\_ref, SourceBioAssays\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- physicalBioAssay\_ref, Target\_assnref-method  
(zzz-DTTCompatibility-class),  
14

physicalBioAssay\_ref,TopLevelBioAssays\_assnreflist-method  
 (zzz-DTTCompatibility-class), 14  
 PhysicalBioAssay\_ref-class  
 (zzz-DTTCompatibility-class), 14  
 physicalBioAssayData\_assnlist  
 (zzz-DTTCompatibility-class), 14  
 physicalBioAssayData\_assnlist,PhysicalBioAssay-method  
 (zzz-DTTCompatibility-class), 14  
 PhysicalBioAssayData\_assnlist-class  
 (zzz-DTTCompatibility-class), 14  
 physicalBioAssaySource\_assnref  
 (zzz-DTTCompatibility-class), 14  
 physicalBioAssaySource\_assnref,FeatureExtraction-method  
 (zzz-DTTCompatibility-class), 14  
 PhysicalBioAssaySource\_assnref-class  
 (zzz-DTTCompatibility-class), 14  
 PhysicalBioAssayTarget\_assnref-class  
 (zzz-DTTCompatibility-class), 14  
 polymerType\_assn  
 (zzz-DTTCompatibility-class), 14  
 polymerType\_assn,BioSequence-method  
 (zzz-DTTCompatibility-class), 14  
 PolymerType\_assn-class  
 (zzz-DTTCompatibility-class), 14  
 position(zzz-DTTCompatibility-class), 14  
 position,Position\_assn-method  
 (zzz-DTTCompatibility-class), 14  
 Position-class  
 (zzz-DTTCompatibility-class), 14  
 position\_assn  
 (zzz-DTTCompatibility-class), 14  
 position\_assn,Feature-method  
 (zzz-DTTCompatibility-class), 14  
 position\_assn,Fiducial-method  
 (zzz-DTTCompatibility-class), 14  
 Position\_assn-class  
 (zzz-DTTCompatibility-class), 14  
 positionDelta  
 (zzz-DTTCompatibility-class), 14  
 positionDelta,PositionDelta\_assn-method  
 (zzz-DTTCompatibility-class), 14  
 PositionDelta-class  
 (zzz-DTTCompatibility-class), 14  
 positionDelta\_assn  
 (zzz-DTTCompatibility-class), 14  
 positionDelta\_assn,FeatureDefect-method  
 (zzz-DTTCompatibility-class), 14  
 positionDelta\_assn,ZoneDefect-method  
 (zzz-DTTCompatibility-class), 14  
 PositionDelta\_assn-class  
 (zzz-DTTCompatibility-class), 14  
 presentAbsent  
 (zzz-DTTCompatibility-class), 14  
 presentAbsent,QuantitationType\_assnlist-method  
 (zzz-DTTCompatibility-class), 14  
 PresentAbsent-class  
 (zzz-DTTCompatibility-class), 14  
 presentAbsent\_ref  
 (zzz-DTTCompatibility-class), 14  
 presentAbsent\_ref,QuantitationType\_assnref-method  
 (zzz-DTTCompatibility-class), 14  
 presentAbsent\_ref,QuantitationTypes\_assnreflist-method  
 (zzz-DTTCompatibility-class), 14  
 presentAbsent\_ref,SourcesQuantitationType\_assnreflist-method  
 (zzz-DTTCompatibility-class), 14



- (zzz-DTTCompatibility-class),  
14
- presentAbsent\_ref,TargetQuantitationType\_assnpropertySets\_assnlist,BioAssayMap-method  
(zzz-DTTCompatibility-class),  
14
- PresentAbsent\_ref-class  
(zzz-DTTCompatibility-class),  
14
- producerTransformation\_assn  
(zzz-DTTCompatibility-class),  
14
- producerTransformation\_assn,DerivedBioAssayDatapropertySets\_assnlist,BioDataCube-method  
(zzz-DTTCompatibility-class),  
14
- ProducerTransformation\_assn-class  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Array-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ArrayManufacture-methodpropertySets\_assnlist,Channel-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ArrayManufactureDeviationpropertySets\_assnlist,CompositeCompositeMap-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Audit-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BibliographicReferencepropertySets\_assnlist,CompositePosition-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BioAssayCreation-methodpropertySets\_assnlist,CompositeSequence-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BioAssayDataCluster-methodpropertySets\_assnlist,CompositeSequenceDimension-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BioAssayDimension-methodpropertySets\_assnlist,Compound-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BioAssayMapping-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BioAssayTreatment-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BioDataTuples-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BioMaterialMeasurement-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BioSample-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BioSequence-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,BioSource-method  
(zzz-DTTCompatibility-class),  
14

- (zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,CompoundMeasurement-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ConcentrationUnit-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Database-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,DatabaseEntry-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,DerivedBioAssay-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,DerivedBioAssayData-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,DerivedSignal-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Description-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,DesignElementMapping-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,DistanceUnit-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Error-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ExpectedValue-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Experiment-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ExperimentalFactor-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ExperimentDesign-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ExternalReference-method  
(zzz-DTTCompatibility-class),  
14
- (zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,FactorValue-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Failed-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Feature-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,FeatureDefect-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,FeatureDimension-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,FeatureExtraction-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,FeatureGroup-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,FeatureInformation-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,FeatureLocation-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,FeatureReporterMap-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Fiducial-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Hardware-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,HardwareApplication-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Hybridization-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Image-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ImageAcquisition-method  
(zzz-DTTCompatibility-class),  
14

- (zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,LabeledExtract-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ManufactureLIMS-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ManufactureLIMSBiomaterial-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,MassUnit-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,MeasuredBioAssay-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,MeasuredBioAssayData-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,MeasuredSignal-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Measurement-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,MismatchInformation-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,NameValueType-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Node-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,NodeContents-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,NodeValue-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,OntologyEntry-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Organization-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Parameter-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ParameterValue-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Person-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,PhysicalArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,PhysicalBioAssay-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Position-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,PositionDelta-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,PresentAbsent-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Protocol-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ProtocolApplication-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,PValue-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,QuantitationTypeDimension-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,QuantitationTypeMap-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,QuantitationTypeMapping-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,QuantityUnit-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Ratio-method

- (zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Reporter-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ReporterCompositeMap-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ReporterDimension-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ReporterGroup-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ReporterPosition-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Security-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,SecurityGroup-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,SeqFeature-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,SeqFeatureLocation-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,SequencePosition-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Software-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,SoftwareApplication-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,SpecializedQuantitationType-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,TemperatureUnit-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,TimeUnit-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Transformation-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Treatment-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,VolumeUnit-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,Zone-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ZoneDefect-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ZoneGroup-method  
(zzz-DTTCompatibility-class),  
14
- propertySets\_assnlist,ZoneLayout-method  
(zzz-DTTCompatibility-class),  
14
- PropertySets\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- protocol (zzz-DTTCompatibility-class),  
14
- protocol,Protocol\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Protocol-class  
(zzz-DTTCompatibility-class),  
14
- protocol\_assnlist  
(zzz-DTTCompatibility-class),  
14
- protocol\_assnlist,Protocol\_package-method  
(zzz-DTTCompatibility-class),  
14
- Protocol\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- protocol\_assnref  
(zzz-DTTCompatibility-class),  
14
- protocol\_assnref,ProtocolApplication-method  
(zzz-DTTCompatibility-class),  
14
- Protocol\_assnref-class  
(zzz-DTTCompatibility-class),

- 14  
 protocol\_package  
 (zzz-DTTCompatibility-class),  
 14
- protocol\_package,MAGE-ML-method  
 (zzz-DTTCompatibility-class),  
 14
- Protocol\_package-class  
 (zzz-DTTCompatibility-class),  
 14
- protocol\_ref  
 (zzz-DTTCompatibility-class),  
 14
- protocol\_ref,Protocol\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14
- Protocol\_ref-class  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplication  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplication,ProtocolApplications\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14
- ProtocolApplication-class  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,ArrayDesign-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,ArrayManufacture-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,BioAssayCreation-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,BioAssayMap-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,BioAssayTreatment-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,CompositeComposition-method  
 (zzz-DTTCompatibility-class),  
 14
- 14  
 protocolApplications\_assnlist,FeatureExtraction-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,FeatureReporterMap-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,Hybridization-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,ImageAcquisition-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,PhysicalArrayDesign-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,QuantitationTypeMap-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,ReporterCompositeMap-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,Transformation-method  
 (zzz-DTTCompatibility-class),  
 14
- protocolApplications\_assnlist,Treatment-method  
 (zzz-DTTCompatibility-class),  
 14
- ProtocolApplications\_assnlist-class  
 (zzz-DTTCompatibility-class),  
 14
- providers\_assnreflist  
 (zzz-DTTCompatibility-class),  
 14
- providers\_assnreflist,Experiment-method  
 (zzz-DTTCompatibility-class),  
 14
- providers\_assnreflist-class  
 (zzz-DTTCompatibility-class),  
 14
- publication  
 (zzz-DTTCompatibility-class),  
 14
- publication,BibliographicReference-method  
 (zzz-DTTCompatibility-class),  
 14
- publication,PublicationList-method  
 (zzz-DTTCompatibility-class),  
 14

- publisher, BibliographicReference-method  
 (zzz-DTTCompatibility-class),  
 14
- pvalue (zzz-DTTCompatibility-class), 14
- pvalue, ConfidenceIndicator\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- pvalue, QuantitationType\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14
- PValue-class  
 (zzz-DTTCompatibility-class),  
 14
- pvalue\_ref  
 (zzz-DTTCompatibility-class),  
 14
- pvalue\_ref, ConfidenceIndicators\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- pvalue\_ref, QuantitationType\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14
- pvalue\_ref, QuantitationTypes\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- pvalue\_ref, SourcesQuantitationType\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14
- pvalue\_ref, TargetQuantitationType\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14
- PValue\_ref-class  
 (zzz-DTTCompatibility-class),  
 14
- quality (zzz-DTTCompatibility-class), 14
- quality, ManufactureLIMS-method  
 (zzz-DTTCompatibility-class),  
 14
- quality, ManufactureLIMSBiomaterial-method  
 (zzz-DTTCompatibility-class),  
 14
- qualityControlDescription\_assn  
 (zzz-DTTCompatibility-class),  
 14
- qualityControlDescription\_assn, ExperimentDesign-method  
 (zzz-DTTCompatibility-class),  
 14
- QualityControlDescription\_assn-class  
 (zzz-DTTCompatibility-class),  
 14
- qualityControlStatistics\_assnlist  
 (zzz-DTTCompatibility-class),  
 14
- qualityControlStatistics\_assnlist, ArrayManufacture-method  
 (zzz-DTTCompatibility-class),  
 14
- qualityControlStatistics\_assnlist, BioSample-method  
 (zzz-DTTCompatibility-class),  
 14
- qualityControlStatistics\_assnlist, BioSource-method  
 (zzz-DTTCompatibility-class),  
 14
- qualityControlStatistics\_assnlist, LabeledExtract-method  
 (zzz-DTTCompatibility-class),  
 14
- QualityControlStatistics\_assnlist-class  
 (zzz-DTTCompatibility-class),  
 14
- quantitationDimension\_assnref  
 (zzz-DTTCompatibility-class),  
 14
- quantitationDimension\_assnref, NodeContents-method  
 (zzz-DTTCompatibility-class),  
 14
- QuantitationDimension\_assnref-class  
 (zzz-DTTCompatibility-class),  
 14
- quantitationType\_assnlist  
 (zzz-DTTCompatibility-class),  
 14
- quantitationType\_assnlist, QuantitationType\_package-method  
 (zzz-DTTCompatibility-class),  
 14
- QuantitationType\_assnlist-class  
 (zzz-DTTCompatibility-class),  
 14
- quantitationType\_assnref  
 (zzz-DTTCompatibility-class),  
 14
- quantitationType\_assnref, QuantitationTypeTuple-method  
 (zzz-DTTCompatibility-class),  
 14
- QuantitationType\_assnref-class  
 (zzz-DTTCompatibility-class),  
 14
- quantitationType\_package

- (zzz-DTTCompatibility-class),  
14
- quantitationType\_package,MAGE-ML-method  
(zzz-DTTCompatibility-class),  
14
- QuantitationType\_package-class  
(zzz-DTTCompatibility-class),  
14
- QuantitationType\_ref-class  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeDimension  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeDimension,QuantitationTypeDimension\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- QuantitationTypeDimension-class  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeDimension\_assnlist  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeDimension\_assnlist,BioAssayData\_package-method  
(zzz-DTTCompatibility-class),  
14
- QuantitationTypeDimension\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeDimension\_assnref  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeDimension\_assnref,DerivedBioAssayData\_package-method  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeDimension\_assnref,MeasuredBioAssayData\_package-method  
(zzz-DTTCompatibility-class),  
14
- QuantitationTypeDimension\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeDimension\_ref  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeDimension\_ref,QuantitationTypeDimension\_ref-assnlist-method  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeDimension\_ref,QuantitationTypeDimension\_ref-assnlist-method  
(zzz-DTTCompatibility-class),  
14
- (zzz-DTTCompatibility-class),  
14
- QuantitationTypeDimension\_ref-class  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeMap  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeMap,QuantitationTypeMap\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- QuantitationTypeMap-class  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeMap\_assnlist,BioAssayData\_package-method  
(zzz-DTTCompatibility-class),  
14
- QuantitationTypeMap\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeMap\_ref  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeMap\_ref,QuantitationTypeMaps\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- QuantitationTypeMap\_ref-class  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeMapping  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeMapping,QuantitationTypeMapping\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- QuantitationTypeMapping-class  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeMapping\_assn  
(zzz-DTTCompatibility-class),  
14
- quantitationTypeMapping\_assn,Transformation-method  
(zzz-DTTCompatibility-class),  
14
- QuantitationTypeMapping\_assn-class  
(zzz-DTTCompatibility-class),  
14

(zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist,DerivedSignal-method  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist,Error-method  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist,ExpectedValue-method  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist,Failed-method  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist,MeasuredSignal-method  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist,PresentAbsent-method  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist,PValue-method  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist,QuantitationTypeMapping-method  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist,Ratio-method  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeMaps\_assnreflist,SpecializedQuantitationType-method  
 (zzz-DTTCompatibility-class),  
 14  
 QuantitationTypeMaps\_assnreflist-class  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypes\_assnreflist  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypes\_assnreflist,QuantitationTypeDimensions-method  
 (zzz-DTTCompatibility-class),  
 14  
 QuantitationTypes\_assnreflist-class  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeTuple  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeTuple,QuantitationTypeTuples\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeTuple-class  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeTuples\_assnlist  
 (zzz-DTTCompatibility-class),  
 14  
 quantitationTypeTuples\_assnlist,DesignElementTuple-method  
 (zzz-DTTCompatibility-class),  
 14  
 QuantitationTypeTuples\_assnlist-class  
 (zzz-DTTCompatibility-class),  
 14  
 quantityUnit  
 (zzz-DTTCompatibility-class),  
 14  
 quantityUnit,Unit\_assn-method  
 (zzz-DTTCompatibility-class),  
 14  
 QuantityUnit-class  
 (zzz-DTTCompatibility-class),  
 14  
 ratio (zzz-DTTCompatibility-class), 14  
 ratio,QuantitationType\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14  
 Ratio-class  
 (zzz-DTTCompatibility-class),  
 14  
 ratio\_ref (zzz-DTTCompatibility-class),  
 14  
 ratio\_ref,QuantitationType\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14  
 ratio\_ref,QuantitationTypes\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14  
 ratio\_ref,QuantitationType\_assnreflist-method  
 (zzz-DTTCompatibility-class),  
 14  
 ratio\_ref,TargetQuantitationType\_assnref-method  
 (zzz-DTTCompatibility-class),  
 14  
 Ratio\_ref-class



- (zzz-DTTCompatibility-class),  
14
- readAnnotation (NetAffxResource-class),  
6
- readAnnotation, NetAffxResource, character, character-method  
(NetAffxResource-class), 6
- readAnnotation, NetAffxResource, character, numeric-method  
(NetAffxResource-class), 6
- readAnnotation, NetAffxResource, missing, AffxAnnotation-class  
(NetAffxResource-class), 6
- readAnnotation, NetAffxResource, numeric, character-method  
(NetAffxResource-class), 6
- readAnnotation, NetAffxResource, numeric, numeric-method  
(NetAffxResource-class), 6
- readArr, 8
- readMage, 9
- readNetAffx, 10
- readXml, 9, 11, 12
- regions\_assnlist  
(zzz-DTTCompatibility-class),  
14
- regions\_assnlist, SeqFeature-method  
(zzz-DTTCompatibility-class),  
14
- Regions\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- releaseDate  
(zzz-DTTCompatibility-class),  
14
- releaseDate, SoftwareApplication-method  
(zzz-DTTCompatibility-class),  
14
- replacedLength  
(zzz-DTTCompatibility-class),  
14
- replacedLength, MismatchInformation-method  
(zzz-DTTCompatibility-class),  
14
- replicateDescription\_assn  
(zzz-DTTCompatibility-class),  
14
- replicateDescription\_assn, ExperimentDesign-method  
(zzz-DTTCompatibility-class),  
14
- ReplicateDescription\_assn-class  
(zzz-DTTCompatibility-class),  
14
- reporter (zzz-DTTCompatibility-class),  
14
- reporter, DesignElement\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- reporter, Reporter\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- reporter\_assnlist  
(zzz-DTTCompatibility-class),  
14
- reporter\_assnlist, DesignElement\_package-method  
(zzz-DTTCompatibility-class),  
14
- Reporter\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- reporter\_assnref  
(zzz-DTTCompatibility-class),  
14
- reporter\_assnref, FeatureReporterMap-method  
(zzz-DTTCompatibility-class),  
14
- reporter\_assnref, ReporterPosition-method  
(zzz-DTTCompatibility-class),  
14
- Reporter\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- reporter\_ref  
(zzz-DTTCompatibility-class),  
14
- reporter\_ref, DesignElement\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- reporter\_ref, Reporter\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- reporter\_ref, Reporters\_assnreflist-method  
(zzz-DTTCompatibility-class),  
14
- Reporter\_ref-class  
(zzz-DTTCompatibility-class),  
14
- reporterCompositeMap

|   |  |
|---|--|
| (zzz-DTTCompatibility-class),<br>14   | (zzz-DTTCompatibility-class),<br>14  |
| reporterCompositeMap,DesignElementMap_assnlist-reported<br>(zzz-DTTCompatibility-class),<br>14        | Dimension_ref<br>(zzz-DTTCompatibility-class),<br>14   |
| reporterCompositeMap,ReporterCompositeMap_assnlist-reported<br>(zzz-DTTCompatibility-class),<br>14    | Dimension_ref,DesignElementDimension_assnref-method<br>(zzz-DTTCompatibility-class),<br>14   |
| ReporterCompositeMap-class<br>(zzz-DTTCompatibility-class),<br>14                                     | ReporterDimension_ref-class<br>(zzz-DTTCompatibility-class),<br>14                           |
| reporterCompositeMap_assnlist<br>(zzz-DTTCompatibility-class),<br>14                                  | reporterGroup<br>(zzz-DTTCompatibility-class),<br>14   |
| reporterCompositeMap_assnlist,DesignElement_package-method<br>(zzz-DTTCompatibility-class),<br>14     | ReporterGroup_assnlist-method<br>(zzz-DTTCompatibility-class),<br>14                         |
| ReporterCompositeMap_assnlist-class<br>(zzz-DTTCompatibility-class),<br>14                            | ReporterGroup-class<br>(zzz-DTTCompatibility-class),<br>14                                   |
| reporterCompositeMap_ref<br>(zzz-DTTCompatibility-class),<br>14                                       | reporterGroup_assnlist<br>(zzz-DTTCompatibility-class),<br>14                                |
| reporterCompositeMap_ref,DesignElementMaps_assnref-list-method<br>(zzz-DTTCompatibility-class),<br>14 | ReporterGroup_assnlist,ArrayDesign_package-method<br>(zzz-DTTCompatibility-class),<br>14     |
| reporterCompositeMap_ref,ReporterCompositeMap_assnref-list<br>(zzz-DTTCompatibility-class),<br>14     | ReporterGroup_assnlist-class<br>(zzz-DTTCompatibility-class),<br>14                          |
| ReporterCompositeMap_ref-class<br>(zzz-DTTCompatibility-class),<br>14                                 | reporterGroup_ref<br>(zzz-DTTCompatibility-class),<br>14                                     |
| reporterCompositeMaps_assnreflist<br>(zzz-DTTCompatibility-class),<br>14                              | reporterGroup_ref,ReporterGroups_assnreflist-method<br>(zzz-DTTCompatibility-class),<br>14   |
| reporterCompositeMaps_assnreflist,CompositeSequences_group<br>(zzz-DTTCompatibility-class),<br>14     | ReporterGroup_ref-class<br>(zzz-DTTCompatibility-class),<br>14                               |
| ReporterCompositeMaps_assnreflist-class<br>(zzz-DTTCompatibility-class),<br>14                        | reporterGroups_assnreflist<br>(zzz-DTTCompatibility-class),<br>14                            |
| reporterDimension<br>(zzz-DTTCompatibility-class),<br>14  | reporterGroups_assnreflist,ArrayDesign-method<br>(zzz-DTTCompatibility-class),<br>14         |
| reporterDimension,DesignElementDimension_assnref-list-method<br>(zzz-DTTCompatibility-class),<br>14   | ReporterGroups_assnreflist,PhysicalArrayDesign-method<br>(zzz-DTTCompatibility-class),<br>14 |
| ReporterDimension-class   | ReporterGroups_assnreflist-class   |

- (zzz-DTTCompatibility-class),  
14
- reporterPosition  
(zzz-DTTCompatibility-class),  
14
- reporterPosition,Coordinate\_assn-method  
(zzz-DTTCompatibility-class),  
14
- reporterPosition,ReporterPositionSources\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ReporterPosition-class  
(zzz-DTTCompatibility-class),  
14
- reporterPositionSources\_assnlist  
(zzz-DTTCompatibility-class),  
14
- reporterPositionSources\_assnlist,ReporterComposition-method  
(zzz-DTTCompatibility-class),  
14
- ReporterPositionSources\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- reporters\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- reporters\_assnreflist,ReporterDimension-method  
(zzz-DTTCompatibility-class),  
14
- reporters\_assnreflist,ReporterGroup-method  
(zzz-DTTCompatibility-class),  
14
- Reporters\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- required(zzz-ARRCompatibility-class),  
13
- required,UserAttribute-method  
(zzz-ARRCompatibility-class),  
13
- roles\_assnlist  
(zzz-DTTCompatibility-class),  
14
- roles\_assnlist,Organization-method  
(zzz-DTTCompatibility-class),  
14
- roles\_assnlist,Person-method  
(zzz-DTTCompatibility-class),  
14
- Roles\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- scale\_assn  
(zzz-DTTCompatibility-class),  
14
- scale\_assn,DerivedSignal-method  
(zzz-DTTCompatibility-class),  
14
- scale\_assn,Error-method  
(zzz-DTTCompatibility-class),  
14
- scale\_assn,ExpectedValue-method  
(zzz-DTTCompatibility-class),  
14
- scale\_assn,Failed-method  
(zzz-DTTCompatibility-class),  
14
- scale\_assn,MeasuredSignal-method  
(zzz-DTTCompatibility-class),  
14
- scale\_assn,NodeValue-method  
(zzz-DTTCompatibility-class),  
14
- scale\_assn,PresentAbsent-method  
(zzz-DTTCompatibility-class),  
14
- scale\_assn,PValue-method  
(zzz-DTTCompatibility-class),  
14
- scale\_assn,Ratio-method  
(zzz-DTTCompatibility-class),  
14
- scale\_assn,SpecializedQuantitationType-method  
(zzz-DTTCompatibility-class),  
14
- Scale\_assn-class  
(zzz-DTTCompatibility-class),  
14
- security(zzz-DTTCompatibility-class),  
14
- security,Security\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Security-class  
(zzz-DTTCompatibility-class),  
14

- [security\\_assnlist](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnlist, AuditAndSecurity\\_package-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [Security\\_assnlist-class](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, Array-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, ArrayDesign-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, ArrayGroup-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, ArrayManufacture-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, Audit-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, BibliographicReference-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, BioAssayCreation-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, BioAssayDataCluster-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, BioAssayDimension-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, BioAssayMap-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, BioAssayTreatment-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, BioSample-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, BioSequence-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, BioSource-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, Channel-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, CompositeCompositeMap-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, CompositeGroup-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, CompositeSequence-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, CompositeSequenceDimension-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, Compound-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, Database-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, DerivedBioAssay-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, DerivedBioAssayData-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, DerivedSignal-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, Description-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, Error-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, ExpectedValue-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref, Experiment-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)

- security\_assnref, ExperimentalFactor-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, ExperimentDesign-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, FactorValue-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Failed-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Feature-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, FeatureDimension-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, FeatureExtraction-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, FeatureGroup-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, FeatureReporterMap-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Fiducial-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Hardware-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, HardwareApplication-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Hybridization-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Image-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, ImageAcquisition-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, LabeledExtract-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, MAGE-ML-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, ManufactureLIMS-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, ManufactureLIMSBiomaterial-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, MeasuredBioAssay-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, MeasuredBioAssayData-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, MeasuredSignal-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Node-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, NodeContents-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Organization-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Parameter-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Person-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, PhysicalArrayDesign-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, PhysicalBioAssay-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, PresentAbsent-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, Protocol-method  
(zzz-DTTCompatibility-class),  
[14](#)
- security\_assnref, ProtocolApplication-method  
(zzz-DTTCompatibility-class),  
[14](#)

- [security\\_assnref,PValue-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,QuantitationTypeDimension-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,QuantitationTypeMap-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,Ratio-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,Reporter-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,ReporterCompositeMap-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,ReporterDimension-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,ReporterGroup-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,Security-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,SecurityGroup-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,SeqFeature-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,Software-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,SoftwareApplication-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,SpecializedQuantitationType-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,Transformation-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,Treatment-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [security\\_assnref,Zone-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [Security-class](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [Security\\_ref](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [Security\\_ref,Security\\_assnref-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [Security\\_ref-class](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroup](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroup,SecurityGroup\\_assnlist-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroup-class](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroup\\_assnlist](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroup\\_assnlist,AuditAndSecurity\\_package-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroup\\_assnlist-class](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroup\\_ref](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroup\\_ref,SecurityGroups\\_assnreflist-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroup\\_ref-class](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroups\\_assnreflist](#)  
 (zzz-DTTCompatibility-class),  
[14](#)
- [SecurityGroups\\_assnreflist,Security-method](#)  
 (zzz-DTTCompatibility-class),  
[14](#)

- SecurityGroups\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- seqFeature  
(zzz-DTTCompatibility-class),  
14
- seqFeature, SeqFeatures\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- SeqFeature-class  
(zzz-DTTCompatibility-class),  
14
- seqFeatureLocation  
(zzz-DTTCompatibility-class),  
14
- seqFeatureLocation, Regions\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- seqFeatureLocation, Subregions\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- SeqFeatureLocation-class  
(zzz-DTTCompatibility-class),  
14
- seqFeatures\_assnlist  
(zzz-DTTCompatibility-class),  
14
- seqFeatures\_assnlist, BioSequence-method  
(zzz-DTTCompatibility-class),  
14
- SeqFeatures\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- sequenceDatabases\_assnlist  
(zzz-DTTCompatibility-class),  
14
- sequenceDatabases\_assnlist, BioSequence-method  
(zzz-DTTCompatibility-class),  
14
- SequenceDatabases\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- sequencePosition  
(zzz-DTTCompatibility-class),  
14
- sequencePosition, Coordinate\_assn-method  
(zzz-DTTCompatibility-class),  
14
- SequencePosition-class  
(zzz-DTTCompatibility-class),  
14
- serialNumber  
(zzz-DTTCompatibility-class),  
14
- serialNumber, HardwareApplication-method  
(zzz-DTTCompatibility-class),  
14
- show, .TypedAssociation-method  
(Internal-class), 4
- show, AffyCompatible-method  
(AffyCompatible-class), 3
- show, NetAffxResource-method  
(NetAffxResource-class), 6
- software (zzz-DTTCompatibility-class),  
14
- software, Software\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Software-class  
(zzz-DTTCompatibility-class),  
14
- software\_assnlist  
(zzz-DTTCompatibility-class),  
14
- software\_assnlist, Protocol\_package-method  
(zzz-DTTCompatibility-class),  
14
- Software\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- software\_assnref  
(zzz-DTTCompatibility-class),  
14
- software\_assnref, SoftwareApplication-method  
(zzz-DTTCompatibility-class),  
14
- Software\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- software\_ref  
(zzz-DTTCompatibility-class),  
14
- software\_ref, Software\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- software\_ref, Softwares\_assnreflist-method

- (zzz-DTTCompatibility-class),  
14
- Software\_ref-class  
(zzz-DTTCompatibility-class),  
14
- softwareApplication  
(zzz-DTTCompatibility-class),  
14
- softwareApplication,SoftwareApplications\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- SoftwareApplication-class  
(zzz-DTTCompatibility-class),  
14
- softwareApplications\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- softwareApplications\_assnreflist,ProtocolApplications\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- SoftwareApplications\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- softwareManufacturers\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- softwareManufacturers\_assnreflist,SoftwareManufacturers\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- SoftwareManufacturers\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- softwares\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- softwares\_assnreflist,Hardware-method  
(zzz-DTTCompatibility-class),  
14
- softwares\_assnreflist,Protocol-method  
(zzz-DTTCompatibility-class),  
14
- softwares\_assnreflist,Software-method  
(zzz-DTTCompatibility-class),  
14
- Softwares\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- sourceBioAssays\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- sourceBioAssays\_assnreflist,BioAssayMap-method  
(zzz-DTTCompatibility-class),  
14
- SourceBioAssays\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- sourceBioMaterialMeasurements\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- sourceBioMaterialMeasurements\_assnreflist,BioAssayCreation-method  
(zzz-DTTCompatibility-class),  
14
- sourceBioMaterialMeasurements\_assnreflist,Hybridization-method  
(zzz-DTTCompatibility-class),  
14
- sourceBioMaterialMeasurements\_assnreflist,Treatment-method  
(zzz-DTTCompatibility-class),  
14
- SourceBioMaterialMeasurements\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- sourceContact\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- sourceContact\_assnreflist,BioSource-method  
(zzz-DTTCompatibility-class),  
14
- SourceContact\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- sourcesQuantitationType\_assnreflist  
(zzz-DTTCompatibility-class),  
14
- sourcesQuantitationType\_assnreflist,QuantitationTypeMap-method  
(zzz-DTTCompatibility-class),  
14
- SourcesQuantitationType\_assnreflist-class  
(zzz-DTTCompatibility-class),  
14
- spacingBetweenCols  
(zzz-DTTCompatibility-class),  
14
- spacingBetweenCols,ZoneLayout-method  
(zzz-DTTCompatibility-class),  
14
- spacingBetweenRows



- (zzz-DTTCompatibility-class),  
14
- spacingBetweenRows, ZoneLayout-method  
(zzz-DTTCompatibility-class),  
14
- spacingsBetweenZonesX  
(zzz-DTTCompatibility-class),  
14
- spacingsBetweenZonesX, ZoneGroup-method  
(zzz-DTTCompatibility-class),  
14
- spacingsBetweenZonesY  
(zzz-DTTCompatibility-class),  
14
- spacingsBetweenZonesY, ZoneGroup-method  
(zzz-DTTCompatibility-class),  
14
- specializedQuantitationType  
(zzz-DTTCompatibility-class),  
14
- specializedQuantitationType, QuantitationType\_ref-method  
(zzz-DTTCompatibility-class),  
14
- SpecializedQuantitationType-class  
(zzz-DTTCompatibility-class),  
14
- specializedQuantitationType\_ref  
(zzz-DTTCompatibility-class),  
14
- specializedQuantitationType\_ref, QuantitationType\_ref-method  
(zzz-DTTCompatibility-class),  
14
- specializedQuantitationType\_ref, QuantitationType\_ref, SeqFeatureLocation-method  
(zzz-DTTCompatibility-class),  
14
- specializedQuantitationType\_ref, SourcesQuantitationType\_ref-method  
(zzz-DTTCompatibility-class),  
14
- specializedQuantitationType\_ref, TargetQuantitationType\_ref-method  
(zzz-DTTCompatibility-class),  
14
- SpecializedQuantitationType\_ref-class  
(zzz-DTTCompatibility-class),  
14
- species\_assn  
(zzz-DTTCompatibility-class),  
14
- species\_assn, BioSequence-method  
(zzz-DTTCompatibility-class),  
14
- species\_assn, CompositeGroup-method  
(zzz-DTTCompatibility-class),  
14
- species\_assn, FeatureGroup-method  
(zzz-DTTCompatibility-class),  
14
- species\_assn, ReporterGroup-method  
(zzz-DTTCompatibility-class),  
14
- Species\_assn-class  
(zzz-DTTCompatibility-class),  
14
- StandardQuantitationType\_ref-class  
(zzz-DTTCompatibility-class),  
14
- startCoord  
(zzz-DTTCompatibility-class),  
14
- startCoord, MatchInformation-method  
(zzz-DTTCompatibility-class),  
14
- strandType  
(zzz-DTTCompatibility-class),  
14
- strandType, SeqFeatureLocation-method  
(zzz-DTTCompatibility-class),  
14
- subregions  
(zzz-DTTCompatibility-class),  
14
- subregions, SeqFeatureLocation-method  
(zzz-DTTCompatibility-class),  
14
- SubregionType\_ref-class  
(zzz-DTTCompatibility-class),  
14
- SubregionType\_ref-method  
(zzz-DTTCompatibility-class),  
14
- substrateType\_assn, ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- SubstrateType\_assn-class  
(zzz-DTTCompatibility-class),  
14
- summaryStatistics\_assnlist

(zzz-DTTCompatibility-class),  
 14  
 summaryStatistics\_assnlist,DerivedBioAssayDataTechnologyType\_assn  
 (zzz-DTTCompatibility-class),  
 14  
 summaryStatistics\_assnlist,MeasuredBioAssayDataTechnologyType\_assn,FeatureGroup-method  
 (zzz-DTTCompatibility-class),  
 14  
 SummaryStatistics\_assnlist-class  
 (zzz-DTTCompatibility-class),  
 14  
 surfaceType\_assn  
 (zzz-DTTCompatibility-class),  
 14  
 surfaceType\_assn,PhysicalArrayDesign-method  
 (zzz-DTTCompatibility-class),  
 14  
 SurfaceType\_assn-class  
 (zzz-DTTCompatibility-class),  
 14  
 target\_assnref  
 (zzz-DTTCompatibility-class),  
 14  
 target\_assnref,BioAssayTreatment-method  
 (zzz-DTTCompatibility-class),  
 14  
 target\_assnref,ImageAcquisition-method  
 (zzz-DTTCompatibility-class),  
 14  
 Target\_assnref-class  
 (zzz-DTTCompatibility-class),  
 14  
 targetQuantitationType\_assnref  
 (zzz-DTTCompatibility-class),  
 14  
 targetQuantitationType\_assnref,Error-method  
 (zzz-DTTCompatibility-class),  
 14  
 targetQuantitationType\_assnref,ExpectedValue-method  
 (zzz-DTTCompatibility-class),  
 14  
 targetQuantitationType\_assnref,PValue-method  
 (zzz-DTTCompatibility-class),  
 14  
 targetQuantitationType\_assnref,QuantitationTypeMap-method  
 (zzz-DTTCompatibility-class),  
 14  
 TargetQuantitationType\_assnref-class  
 (zzz-DTTCompatibility-class),  
 14  
 TechnologyType\_assn  
 (zzz-DTTCompatibility-class),  
 14  
 TechnologyType\_assn-class  
 (zzz-DTTCompatibility-class),  
 14  
 temperatureUnit  
 (zzz-DTTCompatibility-class),  
 14  
 temperatureUnit,Unit\_assn-method  
 (zzz-DTTCompatibility-class),  
 14  
 TemperatureUnit-class  
 (zzz-DTTCompatibility-class),  
 14  
 TemplateFile-class  
 (zzz-ARRCompatibility-class),  
 13  
 timeUnit (zzz-DTTCompatibility-class),  
 14  
 timeUnit,Unit\_assn-method  
 (zzz-DTTCompatibility-class),  
 14  
 TimeUnit-class  
 (zzz-DTTCompatibility-class),  
 14  
 tolerance (zzz-DTTCompatibility-class),  
 14  
 tolerance,ArrayManufacture-method  
 (zzz-DTTCompatibility-class),  
 14  
 tollFreePhone  
 (zzz-DTTCompatibility-class),  
 14  
 tollFreePhone,Organization-method  
 (zzz-DTTCompatibility-class),  
 14  
 tollFreePhone,Person-method  
 (zzz-DTTCompatibility-class),  
 14  
 topLevelBioAssays\_assnreflist  
 (zzz-DTTCompatibility-class),  
 14

topLevelBioAssays\_assnreflist, ExperimentDesign-method  
 (zzz-DTTCompatibility-class),  
 14

TopLevelBioAssays\_assnreflist-class  
 (zzz-DTTCompatibility-class),  
 14

transformation  
 (zzz-DTTCompatibility-class),  
 14

transformation, ProducerTransformation\_assn-method  
 (zzz-DTTCompatibility-class),  
 14

Transformation-class  
 (zzz-DTTCompatibility-class),  
 14

treatment (zzz-DTTCompatibility-class),  
 14

treatment, Treatments\_assnlist-method  
 (zzz-DTTCompatibility-class),  
 14

Treatment-class  
 (zzz-DTTCompatibility-class),  
 14

treatments\_assnlist  
 (zzz-DTTCompatibility-class),  
 14

treatments\_assnlist, BioSample-method  
 (zzz-DTTCompatibility-class),  
 14

treatments\_assnlist, BioSource-method  
 (zzz-DTTCompatibility-class),  
 14

treatments\_assnlist, LabeledExtract-method  
 (zzz-DTTCompatibility-class),  
 14

Treatments\_assnlist-class  
 (zzz-DTTCompatibility-class),  
 14

type (zzz-DTTCompatibility-class), 14

type, ArraySetFile-method  
 (zzz-ARRCompatibility-class),  
 13

type, Measurement-method  
 (zzz-DTTCompatibility-class),  
 14

type, NameValueType-method  
 (zzz-DTTCompatibility-class),  
 14

type, PhysicalArray-method  
 (zzz-ARRCompatibility-class),  
 13

type, TemplateFile-method  
 (zzz-ARRCompatibility-class),  
 13

type, UserAttribute-method  
 (zzz-ARRCompatibility-class),  
 13

type\_assn (zzz-DTTCompatibility-class),  
 14

type\_assn, BioSample-method  
 (zzz-DTTCompatibility-class),  
 14

type\_assn, BioSequence-method  
 (zzz-DTTCompatibility-class),  
 14

type\_assn, DatabaseEntry-method  
 (zzz-DTTCompatibility-class),  
 14

type\_assn, DerivedBioAssay-method  
 (zzz-DTTCompatibility-class),  
 14

type\_assn, Hardware-method  
 (zzz-DTTCompatibility-class),  
 14

type\_assn, NodeValue-method  
 (zzz-DTTCompatibility-class),  
 14

type\_assn, Protocol-method  
 (zzz-DTTCompatibility-class),  
 14

type\_assn, Software-method  
 (zzz-DTTCompatibility-class),  
 14

Type\_assn-class  
 (zzz-DTTCompatibility-class),  
 14

types\_assnlist  
 (zzz-DTTCompatibility-class),  
 14

types\_assnlist, CompositeGroup-method  
 (zzz-DTTCompatibility-class),  
 14

types\_assnlist, ExperimentDesign-method  
 (zzz-DTTCompatibility-class),  
 14

types\_assnlist, FeatureGroup-method

- (zzz-DTTCompatibility-class),  
14
- types\_assnlist,ReporterGroup-method  
(zzz-DTTCompatibility-class),  
14
- Types\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- unit\_assn (zzz-DTTCompatibility-class),  
14
- unit\_assn,Measurement-method  
(zzz-DTTCompatibility-class),  
14
- Unit\_assn-class  
(zzz-DTTCompatibility-class),  
14
- unitName (zzz-DTTCompatibility-class),  
14
- unitName,ConcentrationUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitName,DistanceUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitName,MassUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitName,QuantityUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitName,TemperatureUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitName,TimeUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitName,VolumeUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitNameCV  
(zzz-DTTCompatibility-class),  
14
- unitNameCV,ConcentrationUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitNameCV,DistanceUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitNameCV,MassUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitNameCV,QuantityUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitNameCV,TemperatureUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitNameCV,TimeUnit-method  
(zzz-DTTCompatibility-class),  
14
- unitNameCV,VolumeUnit-method  
(zzz-DTTCompatibility-class),  
14
- upperLeftX  
(zzz-DTTCompatibility-class),  
14
- upperLeftX,Zone-method  
(zzz-DTTCompatibility-class),  
14
- upperLeftY  
(zzz-DTTCompatibility-class),  
14
- upperLeftY,Zone-method  
(zzz-DTTCompatibility-class),  
14
- uri (zzz-DTTCompatibility-class), 14
- uri,BibliographicReference-method  
(zzz-DTTCompatibility-class),  
14
- uri,Database-method  
(zzz-DTTCompatibility-class),  
14
- uri,DatabaseEntry-method  
(zzz-DTTCompatibility-class),  
14
- uri,Description-method  
(zzz-DTTCompatibility-class),  
14
- uri,Hardware-method  
(zzz-DTTCompatibility-class),  
14
- uri,Image-method  
(zzz-DTTCompatibility-class),  
14
- uri,Organization-method  
(zzz-DTTCompatibility-class),

- 14
- uri, Person-method  
(zzz-DTTCompatibility-class),  
14
- uri, Protocol-method  
(zzz-DTTCompatibility-class),  
14
- uri, Software-method  
(zzz-DTTCompatibility-class),  
14
- user (accessors), 2
- user, NetAffxResource-method  
(NetAffxResource-class), 6
- userAttribute  
(zzz-ARRCompatibility-class),  
13
- userAttribute, UserAttributes-method  
(zzz-ARRCompatibility-class),  
13
- UserAttribute-class  
(zzz-ARRCompatibility-class),  
13
- userAttributes  
(zzz-ARRCompatibility-class),  
13
- userAttributes, ArraySetFile-method  
(zzz-ARRCompatibility-class),  
13
- userAttributes, TemplateFile-method  
(zzz-ARRCompatibility-class),  
13
- UserAttributes-class  
(zzz-ARRCompatibility-class),  
13
- userAttributeValue  
(zzz-ARRCompatibility-class),  
13
- userAttributeValue, UserAttribute-method  
(zzz-ARRCompatibility-class),  
13
- UserAttributeValue-class  
(zzz-ARRCompatibility-class),  
13
- value (zzz-ARRCompatibility-class), 13
- value, Control-method  
(zzz-ARRCompatibility-class),  
13
- value, Datum-method  
(zzz-DTTCompatibility-class),  
14
- value, Measurement-method  
(zzz-DTTCompatibility-class),  
14
- value, NameValueType-method  
(zzz-DTTCompatibility-class),  
14
- value, NodeValue-method  
(zzz-DTTCompatibility-class),  
14
- value, OntologyEntry-method  
(zzz-DTTCompatibility-class),  
14
- value, ParameterValue-method  
(zzz-DTTCompatibility-class),  
14
- value\_assn  
(zzz-DTTCompatibility-class),  
14
- value\_assn, FactorValue-method  
(zzz-DTTCompatibility-class),  
14
- Value\_assn-class  
(zzz-DTTCompatibility-class),  
14
- vector, 4
- version (zzz-ARRCompatibility-class), 13
- version, ArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- version, ArraySetFile-method  
(zzz-ARRCompatibility-class),  
13
- version, Database-method  
(zzz-DTTCompatibility-class),  
14
- version, PhysicalArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- version, SoftwareApplication-method  
(zzz-DTTCompatibility-class),  
14
- version, TemplateFile-method  
(zzz-ARRCompatibility-class),  
13
- volume (zzz-DTTCompatibility-class), 14
- volume, BibliographicReference-method

- (zzz-DTTCompatibility-class),  
14
- volumeUnit  
(zzz-DTTCompatibility-class),  
14
- volumeUnit,Unit\_assn-method  
(zzz-DTTCompatibility-class),  
14
- VolumeUnit-class  
(zzz-DTTCompatibility-class),  
14
- warningType\_assn  
(zzz-DTTCompatibility-class),  
14
- warningType\_assn,Reporter-method  
(zzz-DTTCompatibility-class),  
14
- WarningType\_assn-class  
(zzz-DTTCompatibility-class),  
14
- width (zzz-DTTCompatibility-class), 14
- width,ArrayGroup-method  
(zzz-DTTCompatibility-class),  
14
- x (zzz-DTTCompatibility-class), 14
- x,Position-method  
(zzz-DTTCompatibility-class),  
14
- xclass, 9, 11, 12
- xmlValue.XMLAttributeValue, 13
- y (zzz-DTTCompatibility-class), 14
- y,Position-method  
(zzz-DTTCompatibility-class),  
14
- year (zzz-DTTCompatibility-class), 14
- year,BibliographicReference-method  
(zzz-DTTCompatibility-class),  
14
- zone (zzz-DTTCompatibility-class), 14
- zone,ZoneLocations\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- Zone-class  
(zzz-DTTCompatibility-class),  
14
- zone\_assnref  
(zzz-DTTCompatibility-class),  
14
- zone\_assnref,Feature-method  
(zzz-DTTCompatibility-class),  
14
- zone\_assnref,ZoneDefect-method  
(zzz-DTTCompatibility-class),  
14
- Zone\_assnref-class  
(zzz-DTTCompatibility-class),  
14
- zone\_ref (zzz-DTTCompatibility-class),  
14
- zone\_ref,Zone\_assnref-method  
(zzz-DTTCompatibility-class),  
14
- Zone\_ref-class  
(zzz-DTTCompatibility-class),  
14
- zoneDefect  
(zzz-DTTCompatibility-class),  
14
- zoneDefect,Adjustments\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ZoneDefect-class  
(zzz-DTTCompatibility-class),  
14
- zoneGroup (zzz-DTTCompatibility-class),  
14
- zoneGroup,ZoneGroups\_assnlist-method  
(zzz-DTTCompatibility-class),  
14
- ZoneGroup-class  
(zzz-DTTCompatibility-class),  
14
- zoneGroups\_assnlist  
(zzz-DTTCompatibility-class),  
14
- zoneGroups\_assnlist,PhysicalArrayDesign-method  
(zzz-DTTCompatibility-class),  
14
- ZoneGroups\_assnlist-class  
(zzz-DTTCompatibility-class),  
14
- zoneLayout  
(zzz-DTTCompatibility-class),

14  
zoneLayout, ZoneLayout\_assn-method  
(zzz-DTTCompatibility-class),  
14  
ZoneLayout-class  
(zzz-DTTCompatibility-class),  
14  
zoneLayout\_assn  
(zzz-DTTCompatibility-class),  
14  
zoneLayout\_assn, ZoneGroup-method  
(zzz-DTTCompatibility-class),  
14  
ZoneLayout\_assn-class  
(zzz-DTTCompatibility-class),  
14  
zoneLocations\_assnlist  
(zzz-DTTCompatibility-class),  
14  
zoneLocations\_assnlist, ZoneGroup-method  
(zzz-DTTCompatibility-class),  
14  
ZoneLocations\_assnlist-class  
(zzz-DTTCompatibility-class),  
14  
zonesPerX (zzz-DTTCompatibility-class),  
14  
zonesPerX, ZoneGroup-method  
(zzz-DTTCompatibility-class),  
14  
zonesPerY (zzz-DTTCompatibility-class),  
14  
zonesPerY, ZoneGroup-method  
(zzz-DTTCompatibility-class),  
14  
zzz-.TypedAssociation-class  
(Internal-class), 4  
zzz-AffyCompatible-class  
(AffyCompatible-class), 3  
zzz-ARRCompatibility-class, 13  
zzz-DTTCompatibility-class, 14  
zzz-NetAffxCompatibility-class, 16